
LABOROTA 20 control

Betriebsanleitung
Instruction Manual
Mode d'Emploi
Instrucciones
de Empleo
Istruzioni per l'uso

**BEDIENFRONT / FRONT PANEL
 PANNEAU FRONTAL DE COMMANDE / FRONTAL DE MANDO
 PANNELLO FRONTALE DI COMANDO**

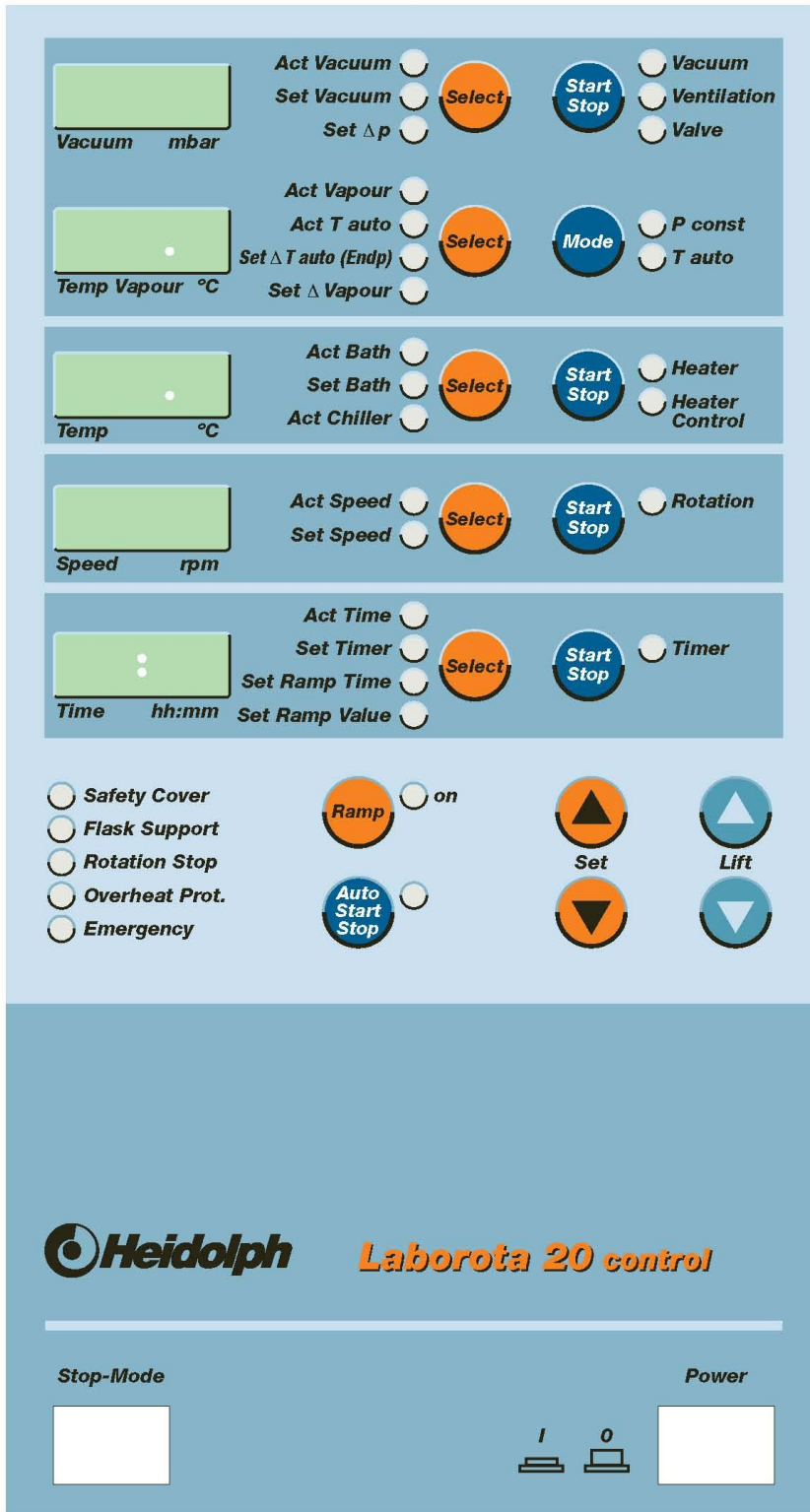DEUTSCH

Seite 4 - 57

ENGLISH

page 58 - 112

FRANCAISE

page 113- 170

ESPAÑOL

página 171 – 227

ITALIANO

página 228 – 284

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EN 61010-1 + EN 61010-2-010

Thank you for buying a Heidolph Instruments product. This unit has been designed, made and inspected in compliance with DIN EN ISO 61010 for long-term and flawless operation.

SUMMARY

BEDIENFRONT / FRONT PANEL	2
SUMMARY	58
PARTS AND ACCESSORIES SUPPLIED	61
GENERAL	62
SAFETY INFORMATION	63
LABOROTA 20 CONTROL, SPECIFICATIONS	64
DISTILLATION SETTING	65
1. General	65
2. Select vacuum control mode	65
3. Solvent data	66
LABORATORY SET-UP	69
SET-UP	70
4. Set-up	70
5. Install vapor tube and PTFE-seal	71
6. Install condensers and expansion vessel	72
7. Install collecting flask in cassette	73
8. Install inlet pipe with valve and screwed cap	74
GLASS-SET -R (CONTROL) 15-300-006-21	76
GLASS-SET -RC (CONTROL) 15-300-006-22	77
GLASS-SET -A (CONTROL) 15-300-006-23	78
GLASS-SET -AC (CONTROL) 15-300-006-24	79
GLASS-SET - A2 (CONTROL) 15-300-006-25	80
GLASS-SET -A2C (CONTROL) 15-300-006-26	81
WORKING WITH THE LABOROTA 20 CONTROL	82
CONTROL PANEL	82
9. Master switch / EMERGENCY-CUT-OUT	82
10. Heating bath lift	83
11. Speed setting	83
12. Heater	84
13. Select vacuum control mode	85
14. p const vacuum control mode	85
15. T auto vacuum control mode	86
16. Calibrate Vacuum pressure sensor	87
17. Auto Start Stop	89

18. Timer.....	90
19. Safety cut-out in an overpressure situation	90
20. Safety cut-off in case of a defect in the vacuum system	92
21. Safety functions	95
TIMER AND RAMP-FUNCTION.....	95
22. Set ramp values	96
23. Display ramp values.....	100
24. Edit ramp values	100
25. Start ramp function.....	100
26. Discontinue ramp function	100
27. Delete ramp	100
INTERFACE	101
28. Connect interface.....	101
29. Interface parameters.....	101
30. RS 232 interface	101
31. RS 232 interface commands.....	101
FUNCTIONS AND HANDLING	102
32. Protective enclosure	102
33. Drive lock	103
34. Tightening of Evaporator Flask	103
35. Flask supports.....	104
36. Automatic water refill (water bath model only)	105
37. Vacuum and water distribution.....	105
38. Pt100 temperature sensor, chiller	106
HEATING BATH FUNCTION AND HANDLING.....	107
39. Fill heating bath (water bath only)	107
40. Drain heating bath (water bath only)	107
41. Condensate drain.....	107
42. Heating bath over temperature protection.....	107
OPTIONS.....	108
43. Vacuum pressure sensor/vacuum valve	108
44. Temperature sensor for vapor.....	108
45. Temperature sensor for T auto vacuum mode.....	108
CLEANING & SERVICING.....	109
UNINSTALL, FORWARD & STORE.....	109
DISPOSAL	109
TROUBLESHOOTING	110
WARRANTY, LIABILITY & COPYRIGHT	111
QUESTIONS / REPAIR WORK.....	111
CE-DECLARATION OF CONFORMITY.....	112

E

Important information

Advise about power cord / mains supply

Caution: mandatory action

Caution: fire- and explosion hazard

Advise about maintenance / repair

PARTS AND ACCESSORIES SUPPLIED

Items contained in your carton (s)

Your LABOROTA 20 carton (s) contain the parts and accessories as listed below. Before disposing of the carton, check the contents for all items on the below parts list.

We recommend to keep packing material for future use.

Item	P/N	qty.	for more details refer to page
LABOROTA 20 control		1	
glass set-R (control)	15-300-006-21	1	page 76
or glass set-RC (control)	15-300-006-22	1	page 77
or glass set-A (control)	15-300-006-23	1	page 78
or glass set-AC (control)	15-300-006-24	1	page 79
or glass set-A2 (control)	15-300-006-25	1	page 80
or glass set-A2C (control)	15-300-006-26	1	page 81
Large coupling	11-300-003-71	1	page 71, (1)
Large split ring	23-30-01-02-31	1	page 71, (13)
PTFE O-Ring	23-30-01-01-11	1	page 71, (5)
PTFE-adapter	23-30-01-02-29-1	1	page 71, (4)
PTFE-seal	23-30-01-01-51-0	1	page 71, (6)
Backing ring	23-30-01-02-28	1	page 71, (7)
Small split ring	23-30-01-02-27	1	page 71, (11)
Small coupling	11-300-003-18	1	page 71, (2)
Vacuum pressure sensor	591-21000-00	1	page 108
Vacuum valve	591-20000-00	1	page 108
Transportation rod	22-30-01-03-17	2	page 70
Cap	11-300-003-28	3	page 70, (5)
Instruction Manual	01-005-004-01	1	

Accessories (optional)

Item	P/N
10 l evaporator flask	15-300-003-16
20 l powder flask	15-300-003-19
10 l powder flask	15-300-003-18
Temperature sensor (for both, boiling temperature and T auto; order 2 ea. when using both options)	14-014-003-06-2
Base cart for LABOROTA 20 control	591-23000-00-1
5 liters bath liquid	515-31000-00

GENERAL

-  Unpack your LR20 carefully. Inspect for damage and report such damage or missing parts to your supplier right away.
-  Read your Instruction Manual carefully. Take time to save time while using your LABOROTA 20. Make sure that every user has read and understood the Instruction Manual.
-  Please keep the Instruction Manual in a place easily accessible to every user.

IF ALL ELSE FAILS, READ THESE INSTRUCTIONS !

For safety reasons, connect your LABOROTA 20 control with a grounded outlet only. (also refer to "Lab Installation")

Using de-ionized or distilled water needs to add 0.2 % of Borax ($\text{Na}_2\text{B}_4\text{O}_7 \cdot 10\text{H}_2\text{O}$) (corrosion inhibitor)

Please note: The use of tapwater as heating media can lead to calcification which can provoke crevice corrosion. Therefore the user is strongly recommended to clean the heating bath with an adequate polish on a regular basis.


SAFETY INFORMATION


- 

Please comply with all safety and accident-prevention regulations, as in force for laboratory work !
- 

Use extra care when working with rotary evaporators.
- 

Use adequate eye protection and protective garments.
- 

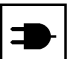Use extra care when working with flammable substances; refer to safety data sheets.
- 
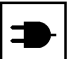
When connecting your unit with your local power supply, please make sure your unit is designed for your local supply voltage; go by data plate on the unit. Please ensure easy access to the power supply.
- 

Turn power switch OFF whenever the unit is not used, or before disconnecting the plug.
- 

Repair work is limited to technicians approved by Heidolph Instruments.
- 

Use extra care when working in the vicinity of flammable and explosive substances. Motors are of non-sparking type, the unit itself however is not explosion-protected.
- 

Caution ! Severe burns may occur when running the heating bath over 50°C. Do not touch hot metal rim of bath, evaporator flask or heating bath fluid.
- 

Never run heating bath dry.
- 

Your LABOROTA 20 needs a solid level stand.
- 


Never attempt to work with damaged or worn-out PTFE gasket (6). Not reaching vacuum pressure wanted may damage your rotary evaporator. Inspect PTFE gasket (6) in regular intervals and replace on condition.
- 

Avoid overpressure in the glassware kit while distilling low-boiling solvents. Aerate glassware kit as required. Overpressure in the glassware kit may damage the evaporator or even blow-up parts of the glassware kit.
- 

In case of disregarding the safety and installation instructions, the safety of the device can be affected.

LABOROTA 20 CONTROL, SPECIFICATIONS

Supply voltage	3 x 400 V 50/60 Hz or 3 x 230 V 50/ 60 Hz 1 x 230V 50/60 Hz
Power rating	4600 VA
Rotation speed (electronic control)	6 - 160 1/min
Heating bath	400 mm dial., stainless steel V2A
Level control, water bath	float switch plus automatic refill valve
Temperature range, water bath	20 - 100 ° C
Temperature range, oil bath	20 - 180 ° C
Heating power	4000 watts
Bath lift	motorized
Over temperature & dry running protection	separate safety circuit
Boil temp. sensor: display range	0 - 150°C
Evaporator flask	20 , 10, or 6 liters optional
Collecting flask	10 liters, features bottom drain valve
Aeration valve	standard
Timer operation	0 – 99:59 h
Setting range, vacuum pressure	1 - 1200 mbar (needs optional vacuum pressure sensor and valve)
Pumping capacity required (vacuum pump)	2.5 - 5 m ³ /h (vacuum pump is optional)
Cooling water consumption	150 - 250 l / h (depending on glass set)
Dimensions (width x depth x height)	970 mm x 620 mm x 1600 mm (with front panel protrusion: 690 mm)
Weight	approx. 120 kg (does not include glass set)
Permissible environment conditions	0-40°C with 80% relative air humidity

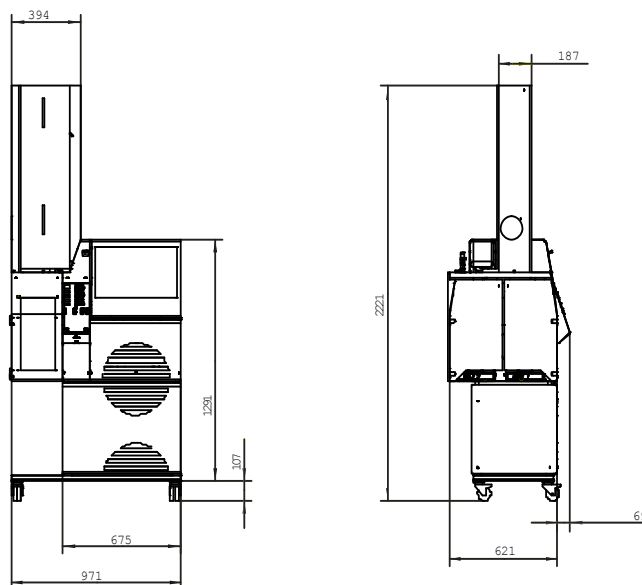DISTILLATION SETTING

1. General

Using rotational evaporators for thermal breakdown of substances needs some basic considerations about optimizing the distillation process. Parameters stipulated below are to be considered for making optimal settings.

- **Rotational speed, evaporator flask**

Increasing rotational speed will increase rate of distillation. Shorter distillation time preserves thermal stability of your substances to be distilled.

- **Difference between heating bath and vapor temperatures:**

This temperature difference should not be less than 20° C to maintain proper rate of distillation. As long as your substances in distillation allow, you may even chose a higher temperature difference.

Rule of thumb: doubling actual temperature difference doubles rate of distillation.

- **Temperature difference between vapor temperature and coolant:**

To maintain adequate condensation rates, we recommend a 20° C temperature difference between the solvent vapor temperature and coolant. A low temperature difference will affect solvent recovery. For substances with high evaporation heat, you should increase temperature difference.

2. Select vacuum control mode

2.1. p_{const} vacuum control mode

The p_{const} mode includes the following features:

- One of the most important modes of control
- Requires optional vacuum valve and vacuum pressure sensor
- Selective distillation of different substances within the scope of their physical properties
- Distillation of solvent mixtures need manual pressure settings
- High reproducibility
- Requires knowledge of thermodynamic substance data

first choice for all standard processes

- For operation & parameter settings refer to page 85
- Pressure (set vacuum) and hysteresis Δp (set Δp) need to be set

E

Select pressure p as to locate boiling point of the substance to be distilled between bath temperature and coolant temperature, make reference to General Notes. For more information refer to table and chart on page 68.

- Δp (hysteresis) value defines max. pressure rise in mbar to open the vacuum valve.

Extended open valve times results in high rates of solvent loss due to extensive gas transfer.

Low-boiling substances require a higher Δp -value (5 – 10 mbar), high-boiling substances a lower value (1 – 5 mbar)

2.2. T_{auto} vacuum control mode

The T_{auto} mode includes the following features:

- Needs optional temperature sensor, vacuum valve and vacuum pressure sensor
- Selective distillation of substances does not need manual pressure setting
- Easy to reproduce
- Easy setting
- No need to know about thermodynamic substance data
-  first choice for low-pollution / zero-loss distillation.
-
- Due to its location, the temperature sensor detects coolant temperature as an approximate only. Set ΔT auto values of 2 – 8°C above coolant temperature for optimum reproducibility. Increasing this temperature span results in a higher rate of distillation with a decrease in reproducibility.

3. Solvent data

- Examples for using table and chart:
- The table displays essential substance data, the chart makes reference to the relationship between pressure and boiling point of some solvents. In this case, $1/T$ is plotted as a function of $\log p$ in accordance with the Clausius-Clapeyron equation.

a) Distillation of a solvent contained in the table:

- on the LABOROTA 20 control set the value from column “vacuum for set point at 40°C” as reference value for Set Vacuum.

b) Distillation of a solvent contained in the chart, with a boiling point other than 40°C:

- Mark the boiling point desired on the temperature axis of the chart
- Draw a line to the right to get an intersection with the solvent
- Vacuum pressure required for this one situation is found by drawing a vertical line from this intersection down to the horizontal axis.

c) Distillation of a solvent not contained in the table or chart:

- The following notes may be helpful to determine the appropriate vacuum pressure:

- The line gradient is a function of enthalpy of evaporation. It is similar for chemically related substances with boiling points close to each other. Hence, the plotted lines may be used as a guide for substances with slightly different boiling points.
- Using a water-jet or diaphragm-type pump, you may lower boiling point by approx. 100°C.
- Rule of thumb: decrease pressure to a half will lower boiling point by approx. 15°C.


Avoid overpressure in the glassware kit while distilling low-boiling solvents. Aerate glassware kit as required. Overpressure in the glassware kit may damage the evaporator or even blow-up parts of the glassware kit.

Solvent	Total formula	MW [g/mol]	MW [°C]	ΔH_{vap} [J/g]	Vacuum for Boiling point at 40°C	
					[mbar]	[mm(Hg)]
Aceton	C ₃ H ₆ O	58,08	56,5	550	556	387
Acetonitril	C ₂ H ₃ N	41,05	81,8	833	230	173
Benzol	C ₆ H ₆	78,11	80,1	549	236	177
n-Butanol (Butylalkohol)	C ₄ H ₁₀ O	74,12	117,5	619	25	19
tert.-Butanol (tert.-Butylalkohol)	C ₄ H ₁₀ O	74,12	82,9	588	130	98
2-Butanon (Methylethylketon)	C ₄ H ₈ O	72,11	79,6	473	243	182
tert.-Butylmethylether	C ₅ H ₁₂ O	88,15	55,0			
Chlorbenzol	C ₆ H ₅ Cl	112,60	132,2	375	36	27
Cyclohexan	C ₆ H ₁₂	84,16	80,7	389	235	176
1,2-Dichlorethan	C ₂ H ₄ Cl ₂	98,96	82,4	336	210	158
1,2-Dichlorethylen (cis)	C ₂ H ₂ Cl ₂	96,94	59,0	320	479	134
1,2-Dichlorethylen (trans)	C ₂ H ₂ Cl ₂	96,94	47,8	313	751	563
Dichlormethan (Methylenchlorid)	CH ₂ Cl ₂	84,93	40,7	373	atm.	atm.
Diethylether	C ₄ H ₁₀ O	74,12	34,6	392	atm.	atm.
Diisopropylether	C ₆ H ₁₄ O	102,20	67,5	318	375	281
Dimethylformamid	C ₃ H ₇ NO	73,09	153,0		11	8
1,4-Dioxan	C ₄ H ₈ O ₂	88,11	101,1	406	107	80
Ethanol	C ₂ H ₆ O	46,07	78,4	879	175	131
Ethylacetat	C ₄ H ₈ O ₂	88,11	77,1	394	240	180
Heptan	C ₇ H ₁₆	85,09	98,4	439	120	90
Hexan	C ₆ H ₁₄	86,18	68,7	370	335	251
Methanol	CH ₄ O	32,04	64,7	1225	337	253
3-Methyl-1-Butanol (Isoamylalkohol)	C ₅ H ₁₂ O	88,15	130,6	593	14	11
Pentachlorethan	C ₂ HCl ₅	202,30	160,5	203	13	10
Pentan	C ₅ H ₁₂	72,15	36,1	382	atm.	atm.
n-Pentanol (Amylalkohol)	C ₅ H ₁₂ O	88,15	137,8	593	11	8
1-Propanol (n-Propylalkohol)	C ₃ H ₈ O	60,10	97,8	787	67	50
2-Propanol (Isopropylalkohol)	C ₃ H ₈ O	60,10	82,5	701	137	103
1,1,2,2-Tetrachlorethan	C ₂ H ₂ Cl ₄	167,90	145,9	247	35	26
Tetrachlorethylen	C ₂ Cl ₄	165,80	120,8	233	53	40
Tetrachlormethan (Carbontetrachlorid)	CCl ₄	153,80	76,7	225	271	203
Tetrahydrofuran	C ₄ H ₈ O	72,11	66,0		357	268
Toluol	C ₇ H ₈	92,14	110,6	425	77	58
1,1,1-Trichlorethan	C ₂ H ₃ Cl ₃	133,40	74,1	251	300	225
Trichlorethylen	C ₂ HCl ₃	131,40	86,7	265	183	137
Trichlormethan (Chloroform)	CHCl ₃	119,40	61,3	263	474	356
Wasser	H ₂ O	18,02	100,0	2259	72	54
Xylol (Isomeren-Gemisch)	C ₈ H ₁₀	106,20	137-143	390	25	19

Conversion factor [mbar] to [mm(Hg)]: $[mm(Hg)] \approx \frac{3}{4} [mbar]$

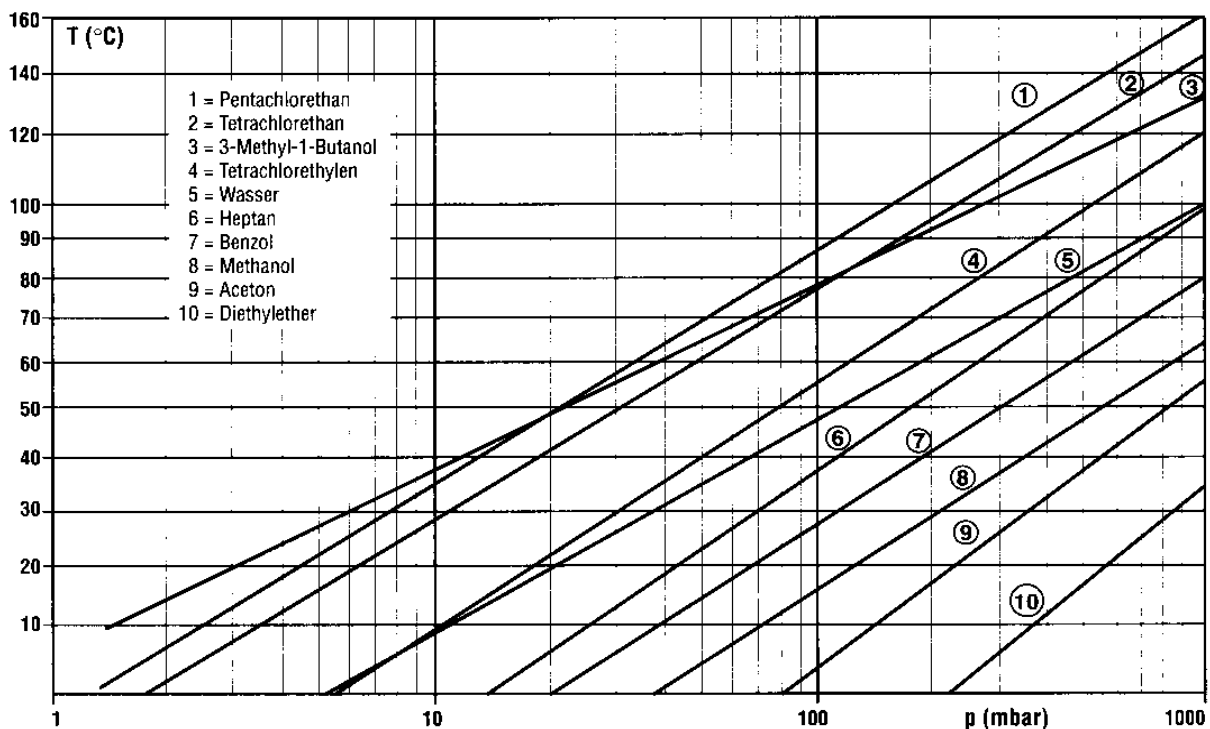LABORATORY SET-UP

SAFETY INFORMATION:

THIS APPLIANCE IS NOT EXPLOSION-PROTECTED. USE EXTREME CARE WHEN USING IN THE VICINITY OF FLAMABLE OR EXPLOSIVE SUBSTANCES.

When connecting this product with your local power supply, make sure you got the right thing ! Check VOLTAGE DATA on DATA PLATE (located in the rear).

This product has to be branched with your local 3-phase network by a skilled electrician.

Wire code:


3 x Phase	GREEN/YELLOW	= protective ground (PE)
	BLUE	= neutral (N)
	BROWN	= phase (L)
	BLACK	= phase (L)
	GREY	= phase (L)

1 x Phase	GREEN	= protective ground (PE)
	WHITE	= neutral (N)
	BLACK	= phase (L)


Carefully unpack all parts, inspect for missing parts by packing list.

4. Set-up

- The **LABOROTA 20 control** with its transportation rods is bolted to the pallet. Remove these screws and carry the unit to its place of installation (**CAUTION: 2 persons needed; in this configuration, unit's weight is about 110 kg**).
- Place the LR 20 on a stable level stand. In ready-to-go configuration, unit's weight will be about 160 kg ! Ideal stand height above ground is 60 cm, this enables easy access to all controls. We recommend the optional **LABOROTA 20 base** (P/N 591-23000-00). Remove transportation rods and keep for later use. Re-install **LABOROTA 20 S control** collecting flask panels (1 & 2) having been removed for transport. Attach panel, to include front door (2) with screws at points (4).
- Through-holes for carrying rods are plugged with caps (5).
- Install protective cabinet (6) on table (7), door pointing towards operator. Use screws (8) pre-installed in attaching holes.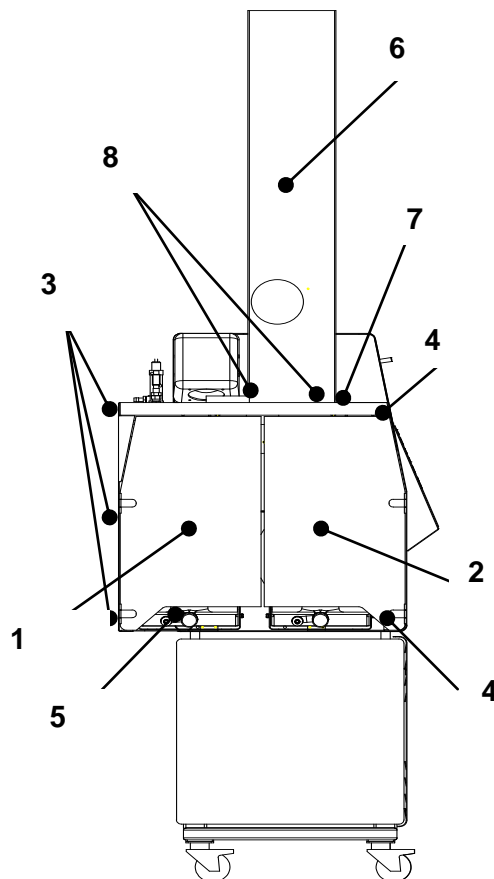

5. Install vapor tube and PTFE-seal

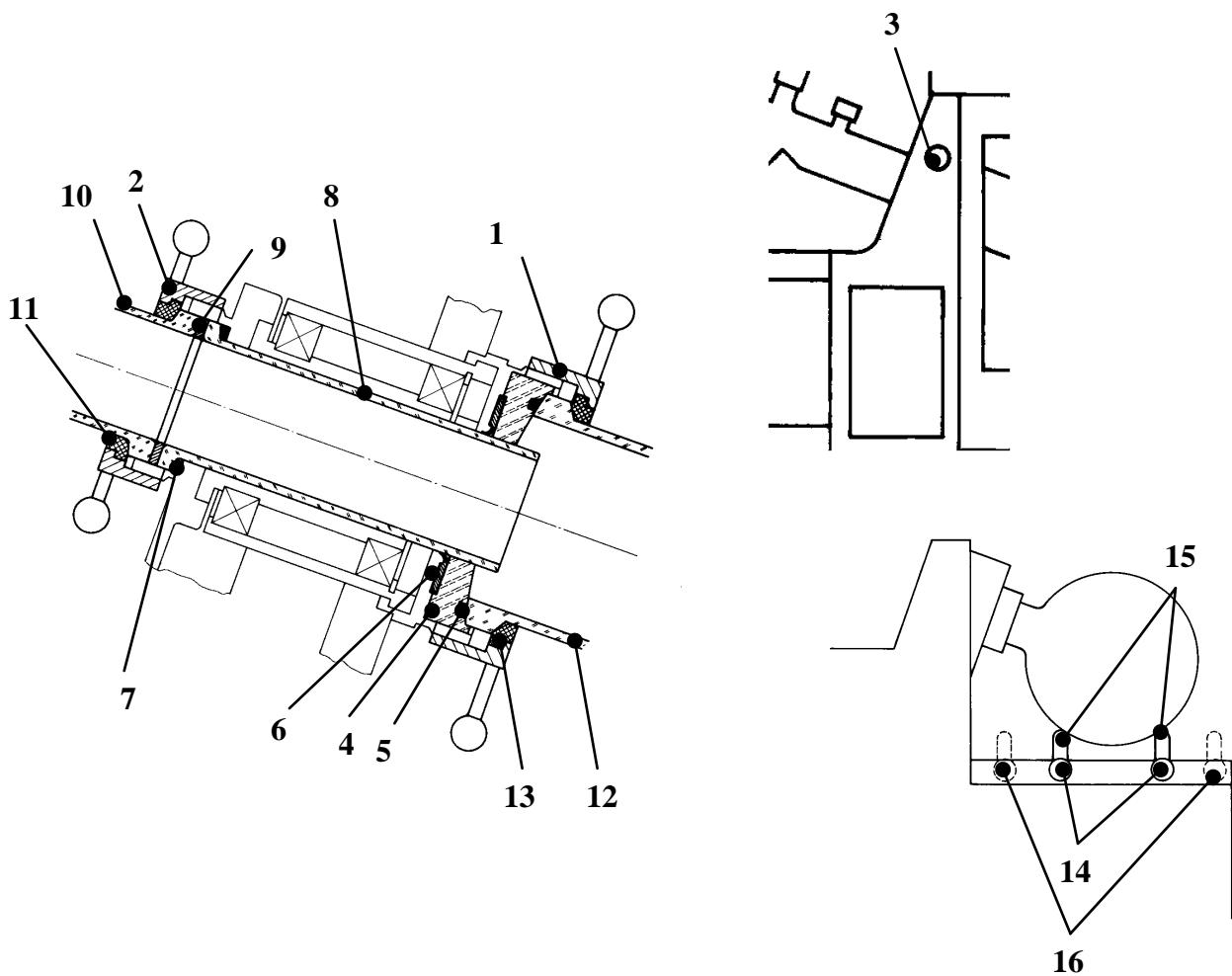Unscrew large-size (1) and small (2) coupling rings. Lock driveshaft by turning drive lock (3) CW to ease loosening and tightening of large-size ring (1).

- Remove PTFE-adapter (4) with built-in PTFE-O-ring (5) and PTFE-seal (6)
- Install backing ring (7) and vapor tube (8) in drive head. Put sandwich gasket (9) on vapor tube coupling flange
- Slide small ring (2) over coupling flange of distributor (10). Expand small split ring (11) and slide over distributor (10) coupling flange, shoulder pointing forward
- Secure assembly (10, 2, 11) tight with small coupling ring (2).

Slide PTFE seal (6) on vapor tube (8), sealing lip pointing inward

Never attempt to work with damaged or worn-out PTFE gasket (6). Not reaching vacuum pressure wanted may damage your rotary evaporator. Inspect PTFE gasket (6) in regular intervals and replace on condition.

- Slide PTFE-adapter (4) with built-in PTFE-O-ring (5) into drive module.
- Slide large-size coupling (1) over evaporator flask (12) coupling flange. Expand large split ring (13) and slide over evaporator flask (12) coupling flange, shoulder pointing forward. Set flask support tubes to 10 liter or 20 liter mark and lock with knobs (14); (I = 10 l, II = 20 l)

E

- Locate evaporator flask on flask lift support tubes (15), tilt flask neck towards drive head, center connecting flange in drive head, so that connecting flange rests against PTFE-O-ring (5) located in PTFE-adapter (4).


Screw large-size ring (1) tight. When evacuating, double-check for firm seat


- Return flask lift support tubes to initial position (16), CAUTION: basic position uses electronic monitor circuit. Leaving this position will freeze all electric functions of the LABOROTA 20 control. This situation is reported by an LED on the control panel (flask support).

6. Install condensers and expansion vessel

- Loosen thumbscrew (2), lift bracket (1) completely and secure with thumbscrew (2)
- Place rubber side of washer (3) on lower bracket (4), line up holes. Put one more washer (3) with rubber side pointing up on upper neck of condenser or expansion vessel (5)
- Plug lower neck of condenser or expansion vessel (5) through cutout in bracket (4) and washer (3), and tilt to an upright position
- Loosen thumbscrew (2) and slide bracket (1) over upper neck of condenser or expansion vessel. Make vertical adjustment with bracket (4)
- Rotate thumbscrew (2) tight when condenser or expansion vessel are in an exactly upright position.

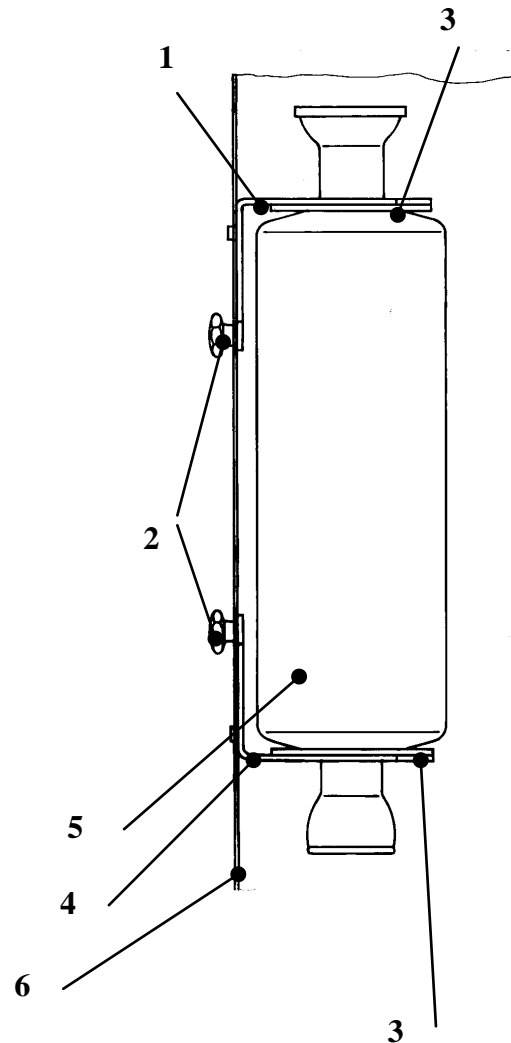
7. Install collecting flask in cassette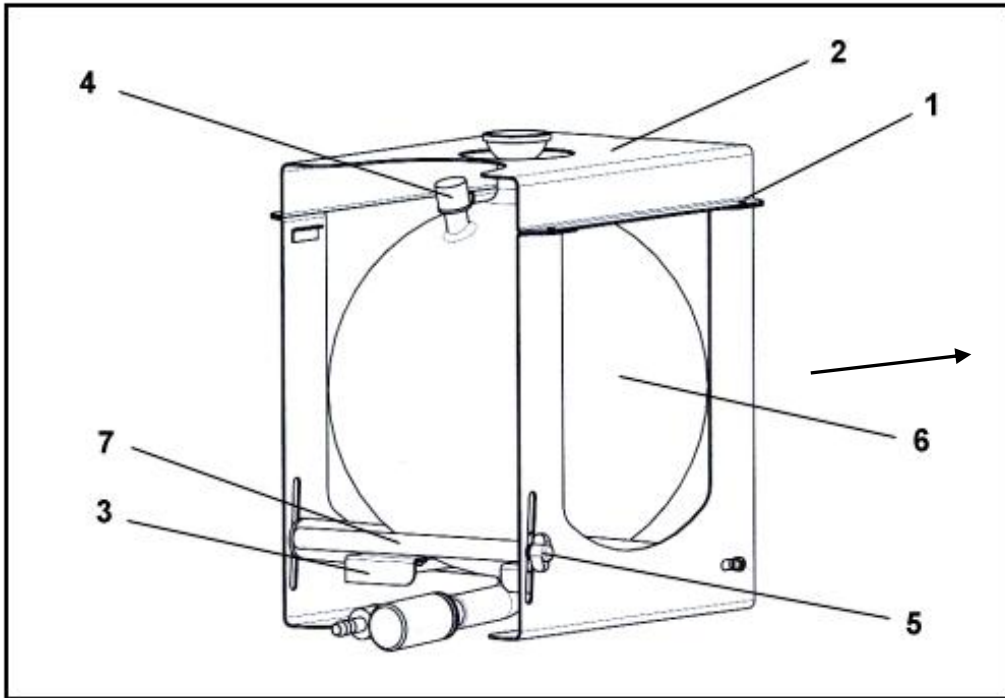


Loosen and remove 4 ea. knurled knobs (1).

- Lift and remove top panel (2).
- Install collecting flask (6) in bottom panel (3). Aerating fitting (4) needs to point forward.
- Reinstall top panel (2) and secure with the 4 knurled knobs (1)
- Install cassette assembly. (with collecting flask) on bracket (8)
- Hold handle (7) of bracket (3), loosen thumbscrew (5), and lower bracket completely; install cassette completely (direction as marked).
- Lift handle (7) until ground joint of collecting flask reaches mating end of condenser, secure with thumbscrew (5)
- use KS45 flask clamp to secure both parts in place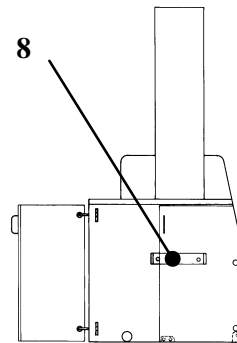


8. Install inlet pipe with valve and screwed cap

1. Valve sealing point is adjusted by turning an Allen screw (4) clockwise with the Allen wrench (3) supplied.
2. Slide threaded ring (5) and gasket kit (5) over PTFE inlet hose (1), as shown. Secure inlet pipe (2) on distributor (6) with threaded ring (5).
3. Install nipple with (7) on inlet pipe (2) with coupling ring (8). Nipple accepts an ID = 6.6 mm hose.

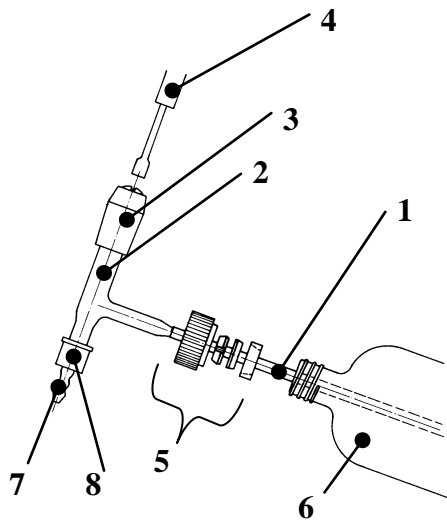
Attention: Never install a valve in main return line of cooling water. A closed valve leads to pressure build-up in the glassware which can lead to damage of glass.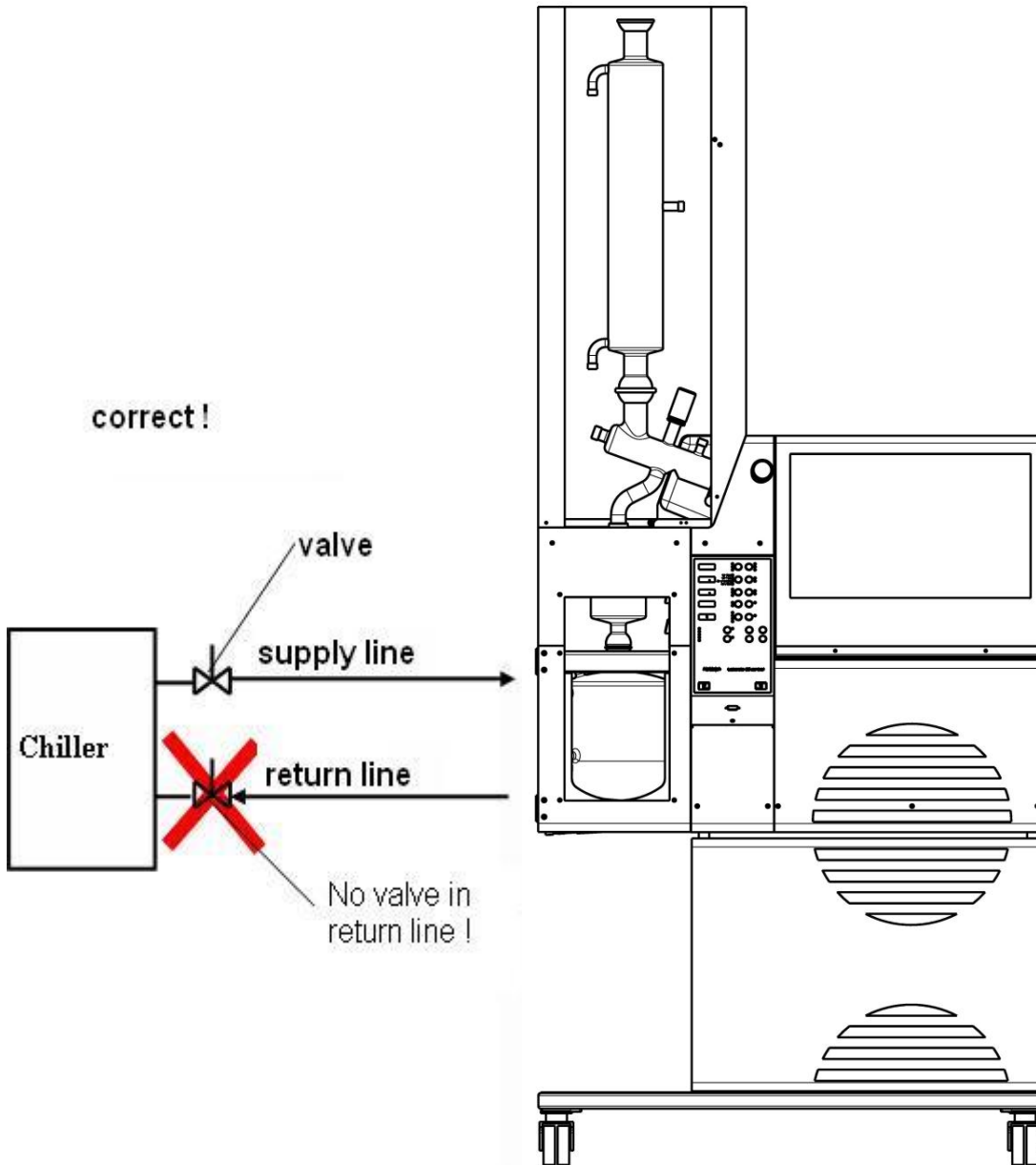

components

A	15-300-001-11	condenser, T auto (R) *
B	15-300-002-31	vacuum fitting
C	15-300-002-58	distributor, includes valve (control)
C1	15-300-002-57	valve GL25/165
E	15-300-002-29	vapor tube, ground
F	15-300-003-17	evaporator flask, 20 l
H	15-300-004-11	collecting flask, 10 l, features bottom drain
H1	15-300-002-56	valve GL25/125 lg.
I	15-300-002-54	inlet pipe
I1	04-03-01-06-02	hose, PTFE 6.5x0.5
N	11-300-003-25	GL25 coupling kit
O	15-300-002-45	valve GL18
P	11-300-005-22	nipple for GL14
Q	23-09-03-01-27	coupling ring GL14
R	23-30-01-02-64	nipple for GL18
S	23-09-03-01-23	coupling ring GL18
T	23-09-03-01-24	screwed cap GL18
U	11-300-002-94	threaded joint KS 64, coated
V	11-300-002-95	flask clamp for S40/S41
W	23-30-01-01-09	sandwich gasket 50 / 69 / 3
X	23-30-01-03-31	spacer sleeve
Z	15-300-001-08	condenser

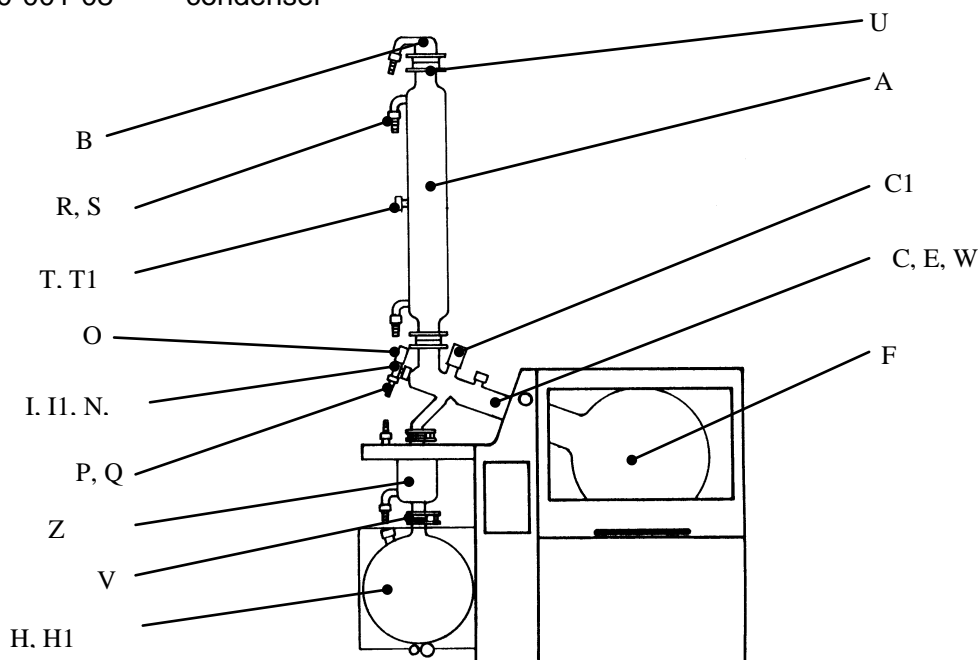
Use this list when checking your shipment for completeness.

* (w. connecting stud for T1 temp. sensor used in T auto vacuum control mode)

GLASS-SET -RC (CONTROL) 15-300-006-22

components

A	15-300-001-11	condenser, T auto (R) *
B	15-300-002-31	vacuum fitting
C	15-300-002-58	distributor, includes valve (control)
C1	15-300-002-57	valve GL25/165
E	15-300-002-29	vapor tube, ground
F	15-300-003-17	evaporator flask, 20 l
G	15-300-002-59	distributor, includes valve
G1	15-300-002-56	valve GL25/125 lg.
H	15-300-004-11	collecting flask, 10 l, features bottom drain
H1	15-300-002-56	valve GL25/125 lg.
I	15-300-002-54	inlet pipe
I1	04-03-01-06-02	hose, PTFE 6.5x0.5
N	11-300-003-25	GL25 coupling kit
O	15-300-002-45	valve GL18
P	11-300-005-22	nipple for GL14
Q	23-09-03-01-27	coupling ring GL14
R	23-30-01-02-64	nipple for GL18
S	23-09-03-01-23	coupling ring GL18
T	23-09-03-01-24	screwed cap GL18
U	11-300-002-94	threaded joint KS 64, coated
V	11-300-002-95	flask clamp for S40/S41
W	23-30-01-01-09	sandwich gasket 50 / 69 / 3
X	23-30-01-03-31	spacer sleeve

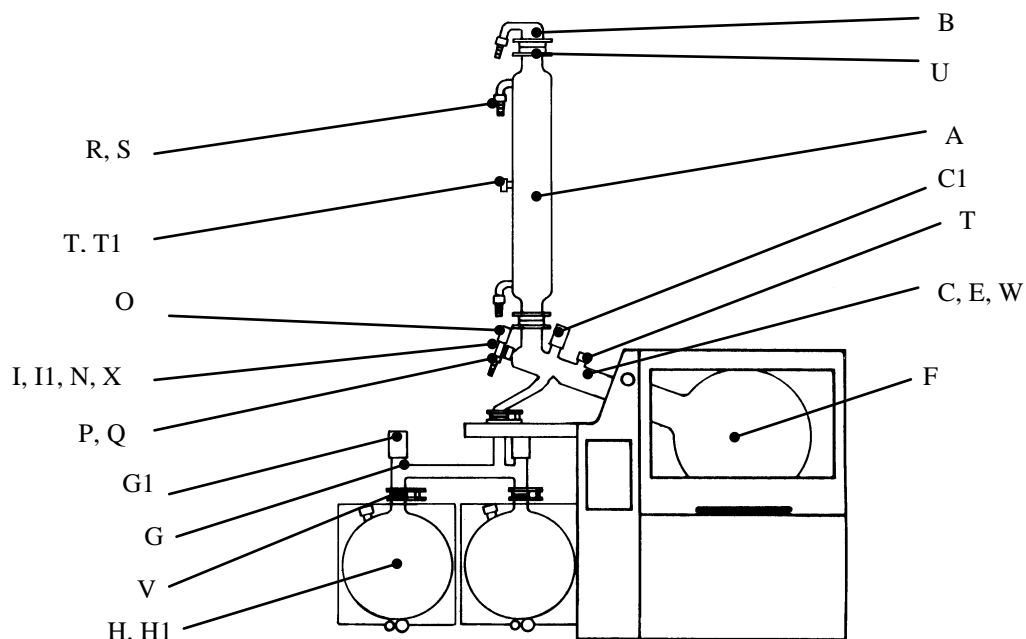
Use this list when checking your shipment for completeness.

* (w. connecting stud for T1 temp. sensor used in T auto vacuum control mode)

GLASS-SET -A (CONTROL) 15-300-006-23

components

A	15-300-001-12	condenser, T auto (A) *
B	15-300-002-31	vacuum fitting
C	15-300-001-09	expansion vessel
D	15-300-002-39	elbow
E	15-300-002-29	vapor tube, ground
F	15-300-003-17	evaporator flask, 20 l
G	15-300-002-40	Tee
H	15-300-004-11	collecting flask, 10 l, features bottom drain
H1	15-300-002-56	valve GL25/125 lg.
I	15-300-002-54	inlet pipe
I1	04-03-01-06-02	hose, PTFE 6.5x0.5
K	15-300-002-41	distributor, to expansion vessel
N	11-300-003-25	GL25 coupling kit
O	15-300-002-45	valve GL18
P	11-300-005-22	nipple for GL14
Q	23-09-03-01-27	coupling ring GL14
R	23-30-01-02-64	nipple for GL18
S	23-09-03-01-23	coupling ring GL18
T	23-09-03-01-24	screwed cap GL18
U	11-300-002-94	threaded joint KS 64, coated
V	11-300-002-95	flask clamp for S40/S41
W	23-30-01-01-09	sandwich gasket 50 / 69 / 3
X	23-30-01-03-31	spacer sleeve
Y	15-300-002-60	valve body & valve
Y1	15-300-002-56	valve GL25 / 125 lg.

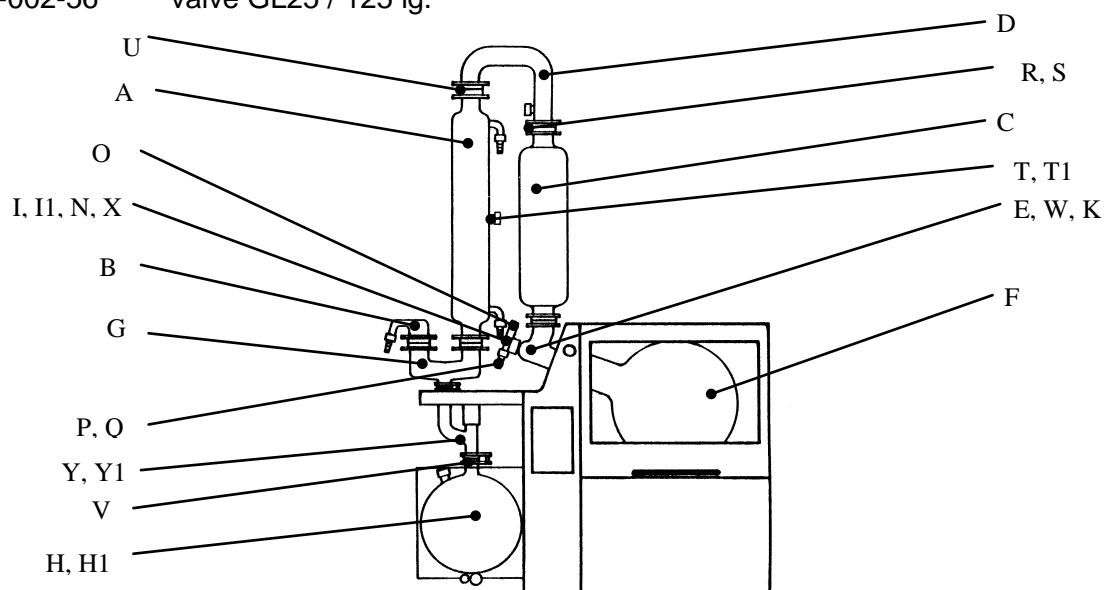
Use this list when checking your shipment for completeness.

* (w. connecting stud for T1 temp. sensor used in T auto vacuum control mode)

GLASS-SET -AC (CONTROL) 15-300-006-24

components

A	15-300-001-12	condenser, T auto (A) *
B	15-300-002-31	vacuum fitting
C	15-300-001-09	expansion vessel
D	15-300-002-39	elbow
E	15-300-002-29	vapor tube, ground
F	15-300-003-17	evaporator flask, 20 l
G	15-300-002-40	Tee
H	15-300-004-11	collecting flask, 10 l, features bottom drain
H1	15-300-002-56	valve GL25/125 lg.
I	15-300-002-54	inlet pipe
I1	04-03-01-06-02	hose, PTFE 6.5x0.5
K	15-300-002-41	distributor, to expansion vessel
L	15-300-002-59	distributor, includes valve
L1	15-300-002-56	valve GL25 / 125lg.
N	11-300-003-25	GL25 coupling kit
O	15-300-002-45	valve GL18
P	11-300-005-22	nipple for GL14
Q	23-09-03-01-27	coupling ring GL14
R	23-30-01-02-64	nipple for GL18
S	23-09-03-01-23	coupling ring GL18
T	23-09-03-01-24	screwed cap GL18
U	11-300-002-94	threaded joint KS 64, coated
V	11-300-002-95	flask clamp for S40/S41
W	23-30-01-01-09	sandwich gasket 50 / 69 / 3
X	23-30-01-03-31	spacer sleeve

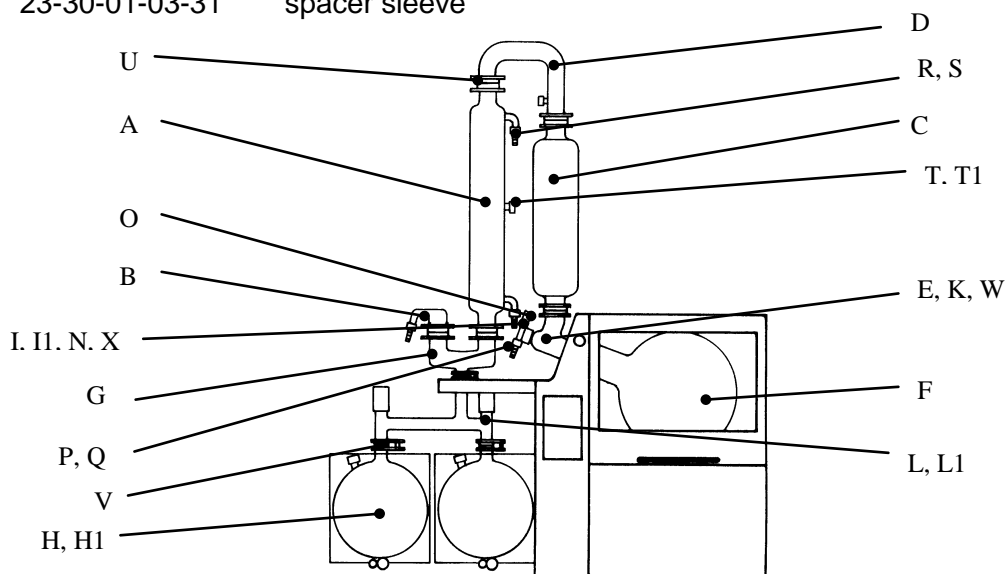
Use this list when checking your shipment for completeness.

* (w. connecting stud for T1 temp. sensor used in T auto vacuum control mode)

GLASS-SET - A2 (CONTROL) 15-300-006-25

components

A	15-300-001-11	condenser, T auto (R) *
A1	15-300-001-12	condenser, T auto (A2) *
B	15-300-002-31	vacuum fitting
C	15-300-001-09	expansion vessel
D	15-300-002-39	elbow
E	15-300-002-29	vapor tube, ground
F	15-300-003-17	evaporator flask, 20 l
G	15-300-002-40	Tee
H	15-300-004-11	collecting flask, 10 l, features bottom drain
H1	15-300-002-56	valve GL25/125 lg.
I	15-300-002-54	inlet pipe
I1	04-03-01-06-02	hose, PTFE 6.5x0.5
K	15-300-002-41	distributor, to expansion vessel
N	11-300-003-25	GL25 coupling kit
O	15-300-002-45	valve GL18
P	11-300-005-22	nipple for GL14
Q	23-09-03-01-27	coupling ring GL14
R	23-30-01-02-64	nipple for GL18
S	23-09-03-01-23	coupling ring GL18
T	23-09-03-01-24	screwed cap GL18
U	11-300-002-94	threaded joint KS 64, coated
V	11-300-002-95	flask clamp for S40/S41
W	23-30-01-01-09	sandwich gasket 50 / 69 / 3
X	23-30-01-03-31	spacer sleeve
Y	15-300-002-60	valve body & valve
Y1	15-300-002-56	valve GL25 / 125 lg.

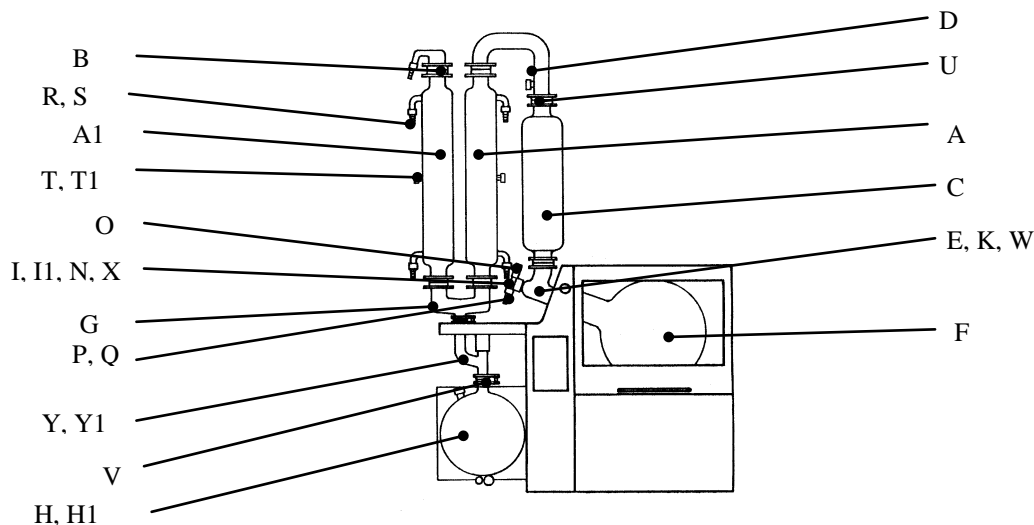
Use this list when checking your shipment for completeness.

* (w. connecting stud for T1 temp. sensor used in T auto vacuum control mode)

GLASS-SET -A2C (CONTROL) 15-300-006-26

components

A	15-300-001-11	condenser, T auto (R) *
A1	15-300-001-12	condenser, T auto (A2) *
B	15-300-002-31	vacuum fitting
C	15-300-001-09	expansion vessel
D	15-300-002-39	elbow
E	15-300-002-29	vapor tube, ground
F	15-300-003-17	evaporator flask, 20 l
G	15-300-002-40	Tee
H	15-300-004-11	collecting flask, 10 l, features bottom drain
H1	15-300-002-56	valve GL25/125 lg.
I	15-300-002-54	inlet pipe
I1	04-03-01-06-02	hose, PTFE 6.5x0.5
K	15-300-002-41	distributor, to expansion vessel
N	11-300-003-25	GL25 coupling kit
O	15-300-002-45	valve GL18
P	11-300-005-22	nipple for GL14
Q	23-09-03-01-27	coupling ring GL14
R	23-30-01-02-64	nipple for GL18
S	23-09-03-01-23	coupling ring GL18
T	23-09-03-01-24	threaded cap GL18
U	11-300-002-94	threaded joint KS 64, coated
V	11-300-002-95	flask clamp for S40/S41
W	23-30-01-01-09	sandwich gasket 50 / 69 / 3
X	23-30-01-03-31	spacer sleeve
Y	15-300-002-59	distributor, includes valve
Y1	15-300-002-56	alve GL25 / 125 lg.

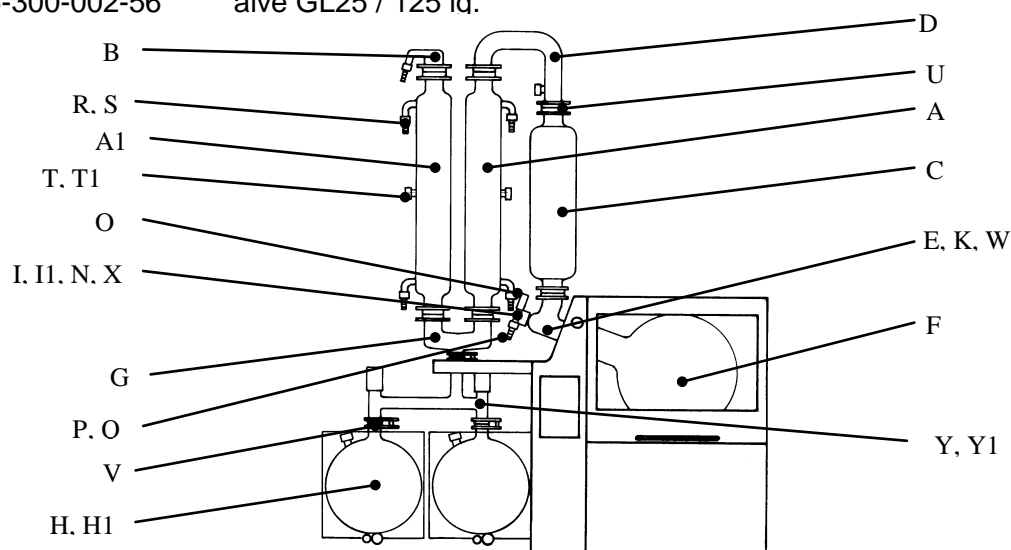Use this list when checking your shipment for completeness.

* (w. connecting stud for T1 temp. sensor used in T auto vacuum control mode)

WORKING WITH THE LABOROTA 20 CONTROL

Use rotary evaporators with appropriate care. Comply with standard laboratory safety practices. Be cautious when working with flammable solvents; motors and limit switches are of non-sparking type but we cannot guarantee that the unit is explosion-proof.

SAFETY INFORMATION:


WHEN CONNECTING YOUR UNIT WITH YOUR LOCAL POWER SUPPLY, PLEASE MAKE SURE YOUR UNIT IS DESIGNED FOR YOUR LOCAL SUPPLY VOLTAGE; GO BY DATA PLATE ON THE UNIT.

CONNECT YOUR ROTARY EVAPORATOR WITH A GROUNDED OUTLET ONLY.

CONTROL PANEL

9. Master switch / EMERGENCY-CUT-OUT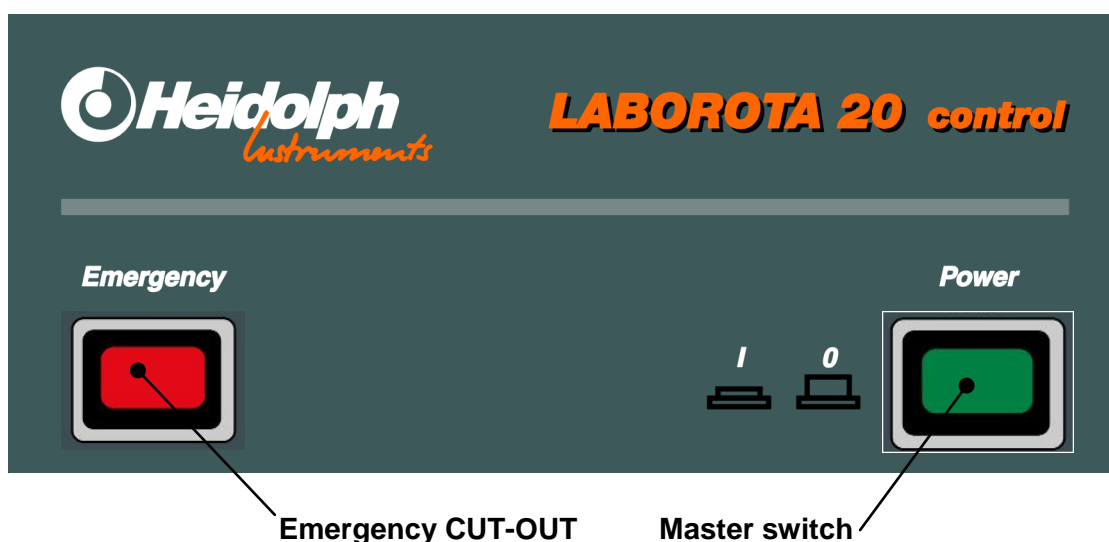

The master switch disconnects all wires from power supply, and serves as a RESET button. (after Emergency CUT-OUT).

The **Emergency CUT-OUT** discontinues operation of the rotary evaporator immediately.

- Stop rotation
- Lowers heating bath
- Opens aerating valve (vacuum valve closes)
- Turns OFF heating circuit
- Buzzer sounds
- Emergency LED turns ON

CAUTION: if entry of ambient air causes unwanted reactions, take adequate action (inert gas, or keep aerating valve closed).

System re-start after an Emergency CUT-OUT situation requires system RESET. Make sure the problem is fixed before resetting. See an expert, if required.

To Reset, turn **Master switch** OFF & ON.

10. Heating bath lift


Heating bath lift RAISES, as long as the **Lift UP** key is held depressed.

Heating bath lift LOWERS, as long as the **Lift DOWN** key is held depressed.

At upper and lower limits, lift motion stops automatically.

The LABOROTA 20 control WB (water bath) features a fully automatic water refill system with float switch.

Exceeding the max. fill mark (about 30 mm below tank rim, monitored by float switch), the **Lift UP** motion is discontinued automatically.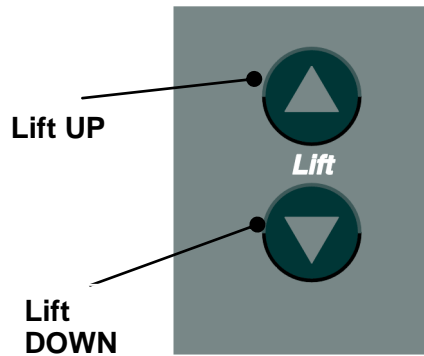


float switch must be connected

As water in the heating bath evaporates, the heating bath itself travels up in small increments for compensation, until reaching the upper stop.

If max fill level is not sensed at the upper stop, the water will be refilled automatically.

The LABOROTA 20 control OB (oil bath) features neither refill or float switch.

11. Speed setting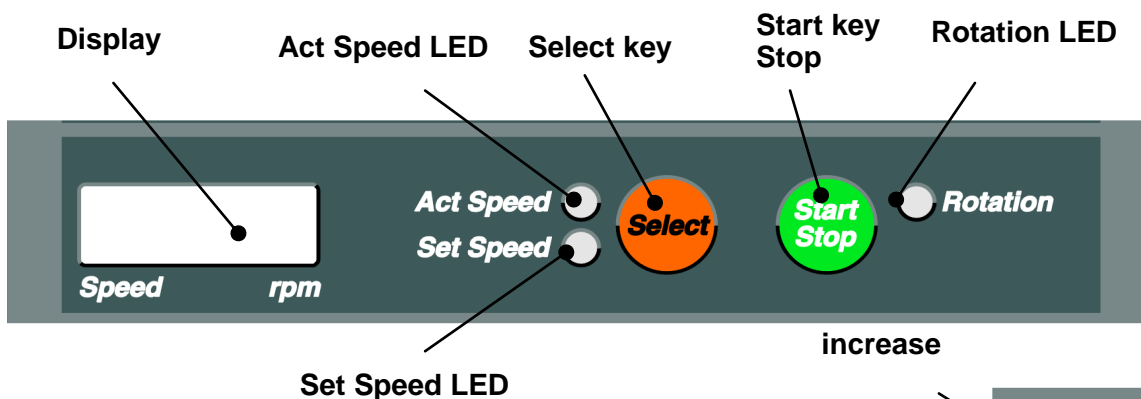

Change between Act Speed and Set Speed display with Select key. **Act Speed LED** or **Set Speed LED** illuminate.

Select **LED Set Speed** for speed setting.

Nominal speed (rpm) appears in the display.

Change nominal speed with **Set keys**.

5 seconds after last entry, display will change to "Act", last entry will be saved.

Start rotation with **Start Stop key**.

Hit this key once more to stop rotation.


CAUTION: for safety reasons, start rotation at low speeds.

Rotational speed can be increased during operation.

Never exceed a speed of 100 rpm, while working with a 20 l evaporation flask.


If rotation does not start with the START/STOP key, please check for illumination of one or more of the Safety Cover, Flask Support or Rotation Stop LED's.

If this occurs check if the enclosure is open, flask support not in one of its extreme positions, or the drive drive lock engaged.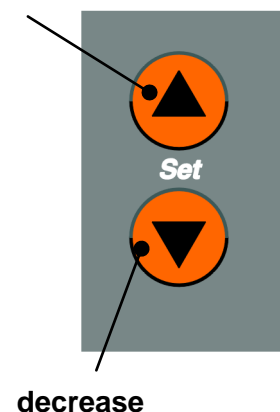

For more information refer to Troubleshooting Section.

12. Heater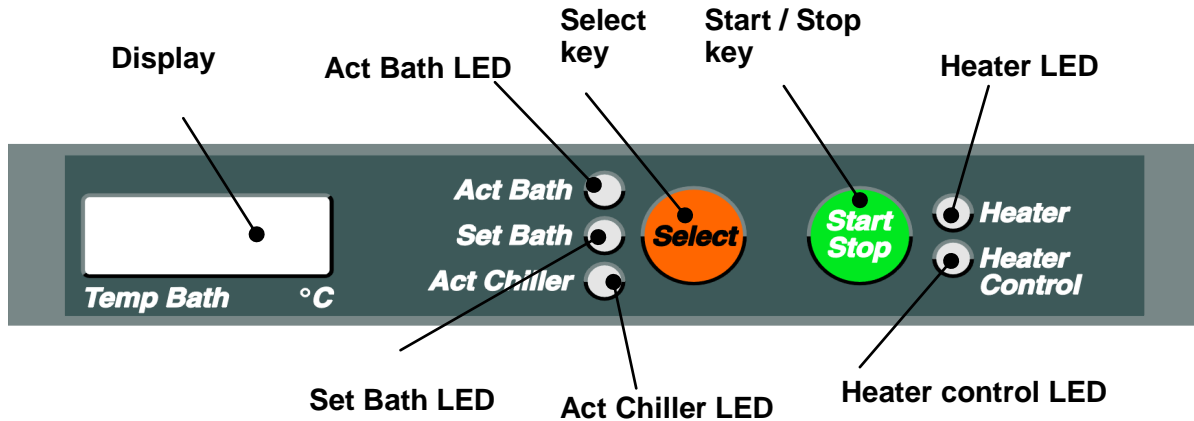

Change between Act Bath, Set Bath and Act Chiller with set key. Act Bath, Set Bath or Act Chiller LED turns ON.

Act Chiller is the actual coolant temperature, which is the basis for the T auto vacuum control mode.

To set heating bath temperature, select **Set Bath LED**.

Display now shows the heating bath's nominal temperature with a 1/10 of a degree C resolution.

Change heating bath temperature with the **Set keys**.

Not pressing one of these keys for 5 seconds will cause the display returns to "Act" mode. The last figure entered is saved.

The **Start Stop** key turns the heater ON.


Pressing the key once more will discontinue heating.

The Heater LED illuminates as long as the heating mode is kept active, the **Heater control LED** illuminates as long as heating current flows. (Act Bath < Set Bath)

Should the Heater LED remain dim after pressing the Start Stop key, please check to see if the Over temperature Protection LED is ON or flashing.

For more information refer to Troubleshooting Section.

increase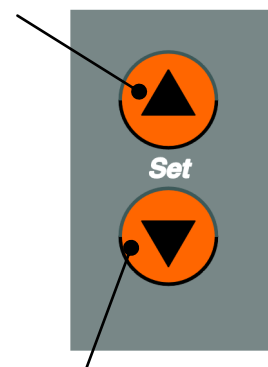

decrease


13. Select vacuum control mode

Vacuum control requires the optional vacuum pressure valve and –sensor.


The **Mode key** changes between p const and T auto vacuum control modes. **p const** or **T auto LED** illuminates.

You may change between these vacuum control modes during operation. For more information about vacuum control modes please refer to Distillation Settings Section.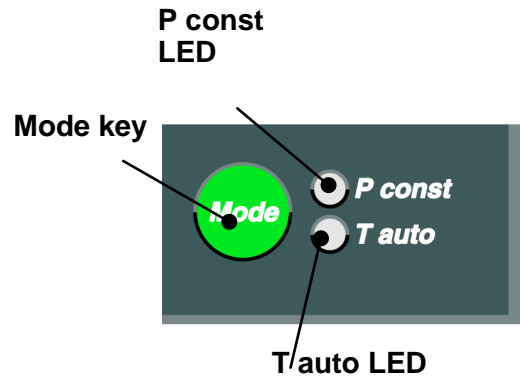

14. p const vacuum control mode

In this mode, a vacuum level once selected will be maintained. **Set Vacuum** key sets the level and **Set Δp** your control span Δp (hysteresis). **Act Vacuum** key provokes display of actual pressure.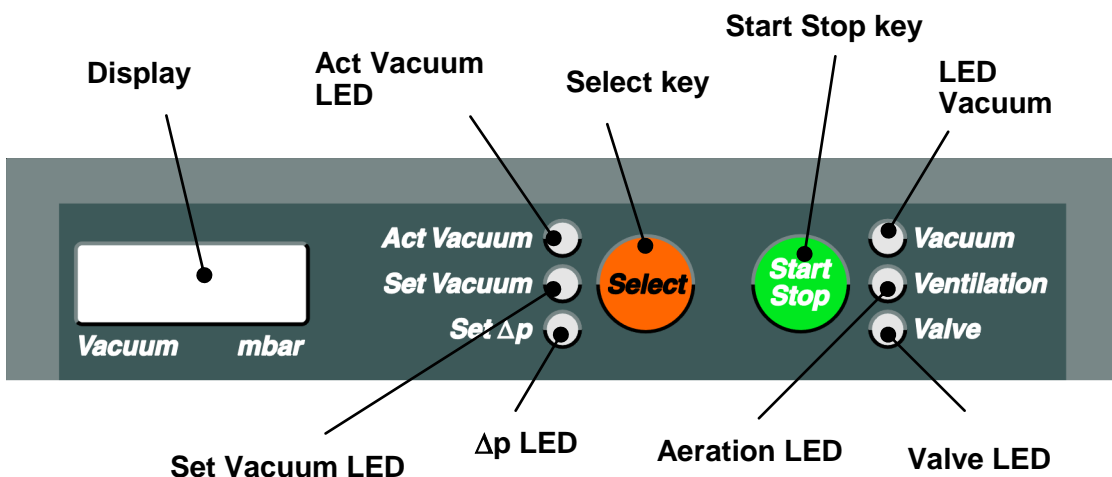

Change between **Act Vacuum**, **Set Vacuum** and **Set Δp** with **Select key**. **Act Vacuum LED**, **Set Vacuum LED**, or **Set Δp LED** will illuminate.

Select **Set Vacuum LED** for setting the nominal vacuum level.

The nominal suction pressure in mbar appears in your **display**.

Change nominal suction pressure level with the **Set keys**.

Not pressing one of these keys for 5 seconds will return the unit to “Act” mode. The last figure entered is saved.


Select Set Δp LED for setting hysteresis of Δp.

Use SET keys as required.

The **Start Stop key** starts evacuation.

Press the key for an instant to interrupt evacuation without opening the aerating valve, Valve LED flashing; one more brief press on this key continues evacuation.

increase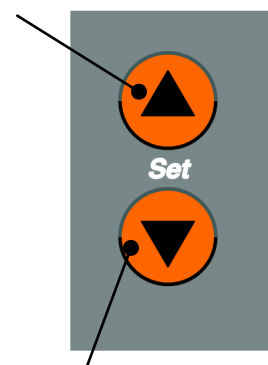

decrease

E

Extended actuation of the Start Stop key discontinues evacuation, the aerating valve opens, the vacuum valve is closed.

Vacuum LED illuminates, when “vacuum control” was selected with the Start Stop key.

Ventilation LED illuminates, when the aerating valve is open.

Valve LED illuminates, when the vacuum valve is open.

Valve LED flashes, when evacuation was interrupted without aeration.

15. T auto vacuum control mode

In this mode, vacuum pressure level is controlled by an optional temperature sensor. For installation refer to Options Section (page 93).

T auto is a low-pollution mode. This special vacuum control allows very low quantities of solvents to be extracted by the vacuum pump.


The T auto temperature sensor detects temperature in the middle of the condenser.

For cooling water temperature, the LABOROTA 20 control uses a separate temperature sensor (act Chiller).

Set ΔT auto is set to 2 – 10 °C (temperature difference from cooling water).

A high Set ΔT auto value results in high-speed distillation, whereas a low value renders a slow and controlled distillation process.

When reaching the temperature difference (in the condenser) once set with Set ΔT auto, the vacuum valve closes; as soon as the temperature drop equals Set ΔV auto (hysteresis), the valve will open once more. A current ΔV auto setting is 0.5°C.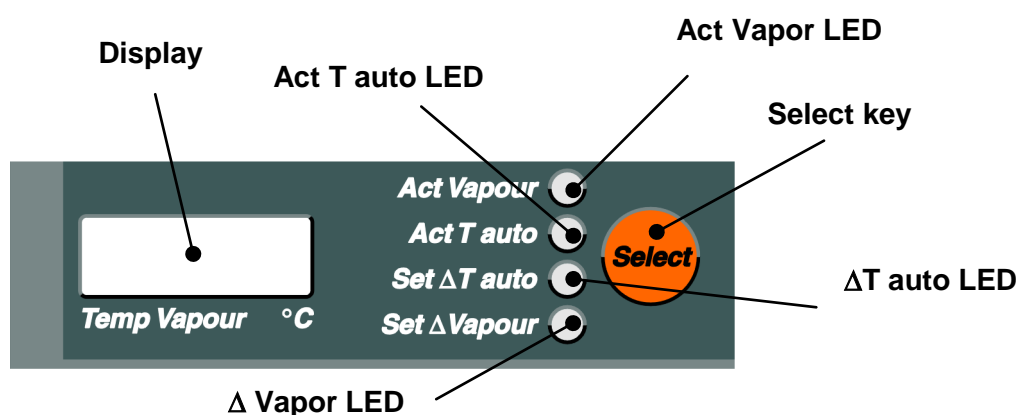


Pressing the Select key changes display reading between **Act Vapor**, **Act T auto**, **Set ΔT auto** and **Set ΔV auto**.

The Act Vapor, Act T auto, Set ΔT auto or Set ΔV auto will illuminate.

Act T auto shows command temperature in the condenser for mode **T auto**, i.e. at point 5 (see page 108), as soon as the temperature sensor is installed and connected with #4 plug connector (T auto temperature sensor) (see page 108).

Act Vapor shows vapor temperature in the

increase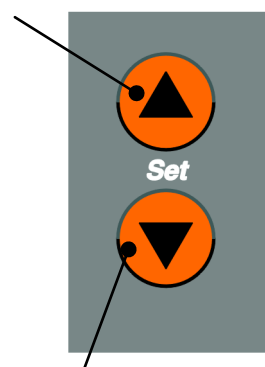


decrease

distributor, i.e. point 1 (see page 93) as soon as the temperature sensor is installed and connected with #3 plug connector (boiling temperature sensor) (see page 93).

For **Set ΔT auto** you need to select the **Set ΔT auto LED**.

Next step is entering final pressure; exceeding this value automatically discontinues distillation. After about 5 seconds, "**Endp**" appears in display II, at this point select final pressure. Final pressure selected appears in the display I.

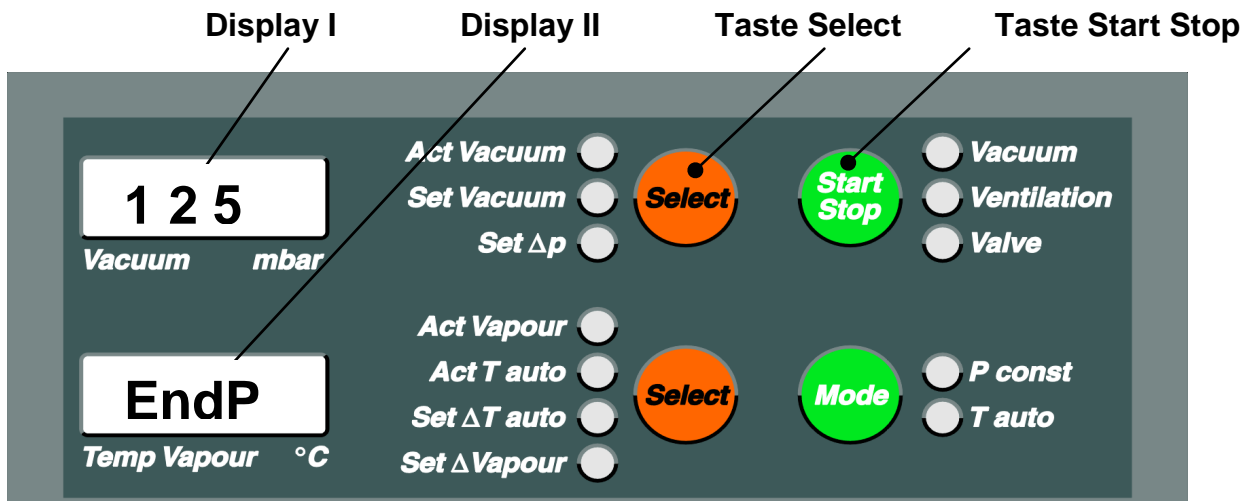Reaching this final pressure during distillation, will discontinue distillation right away. Also refer to Auto Start Stop function (see page 89)


If distillation should not be discontinued, enter "1" in "**Endp** mode", which never can be reached by physical condition.

For hysteresis setting select **LED Δ Vapor**.

Not pressing a key for 5 seconds, display returns to "**Act**" mode. The last figure entered is saved. Start evacuation, as described in page 85.


16. Calibrate Vacuum pressure sensor

- Vacuum pressure sensor and vacuum valve must be installed on the LABOROTA 20 control and a vacuum pump connected.
- A separate, calibrated vacuum gauge should be installed in the vacuum line near condenser
- Turn vacuum pump ON
- Turn LABOROTA 20 control master switch (**Power**) OFF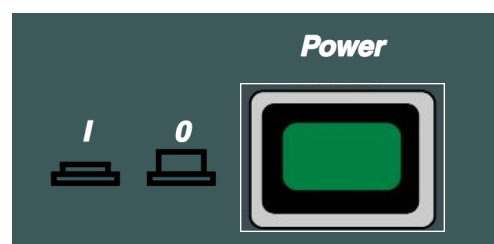


E

- Depress Select and Start Stop keys near Display I together and hold for 5 seconds, turn master switch ON, release keys
- „CAL“ is shown in the **Display Speed**, „VAC“ in the **Display Time**
- 750 is shown in Display I and 1 is shown in Display II (1st step).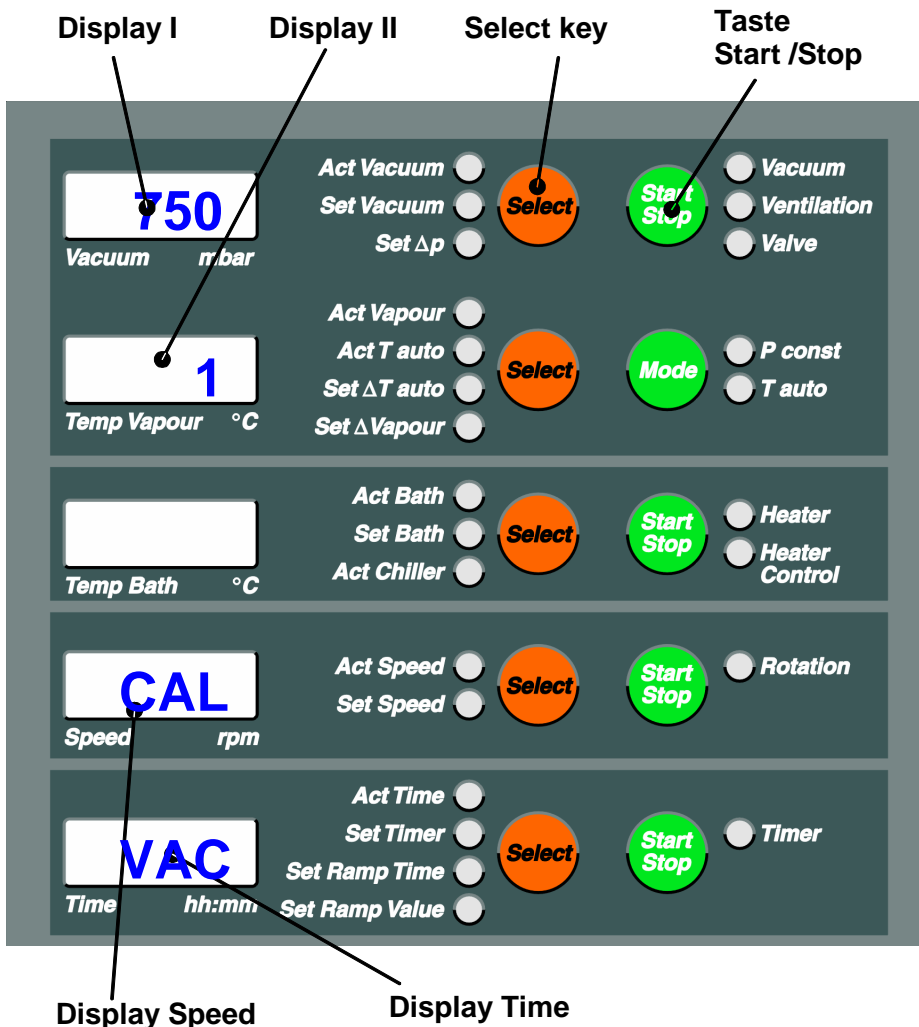

- Now, set the pressure measured with the separate vacuum gauge in Display I, by pressing the Set increase key or Set decrease key
- Confirm with **Start Stop** key located next to Display I
- Display I to show 50 and display II 2 (2nd step).
- At this point, the LABOROTA 20 control automatically opens the vacuum valve and evacuates until reaching vacuum limit (the value on the separate vacuum gauge won't change anymore)
- In Display I, set the pressure measured with the separate vacuum gauge in Display I, by pressing the Set increase key or Set decrease key
- Confirm with **Start Stop** key located next to Display I. As confirmation for the accepted calibration values, „0“ appears in Display I and Display II
- Turn master switch (**Power**) OFF and ON
- Calibration is complete

increase

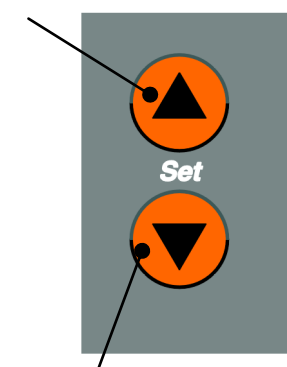decrease

17. Auto Start Stop

The Auto Start Stop key combines several functions; this way, distillation may be started and stopped automatically. Before such action, enter all distillation parameters and start the following functions with the Start Stop key:

- Lift UP
- Rotation ON
- Heater ON
- Vacuum ON
- Start timer (LED Act Time flashing)

An LED beside the Auto Start Stop key confirms such action.

CAUTION: if one or more functions are already active, e.g. heater, only the remaining functions will be started.


The Time display shows time elapsed from pressing the Auto Start Stop key.

Pressing the Auto Start Stop key once more will reset the functions started:

- Lift DOWN
- Rotation OFF
- Heater OFF
- Vacuum OFF
- Buzzer sounds for 5 seconds
-

During this automatic mode, all distillation parameters may be changed!


18. Timer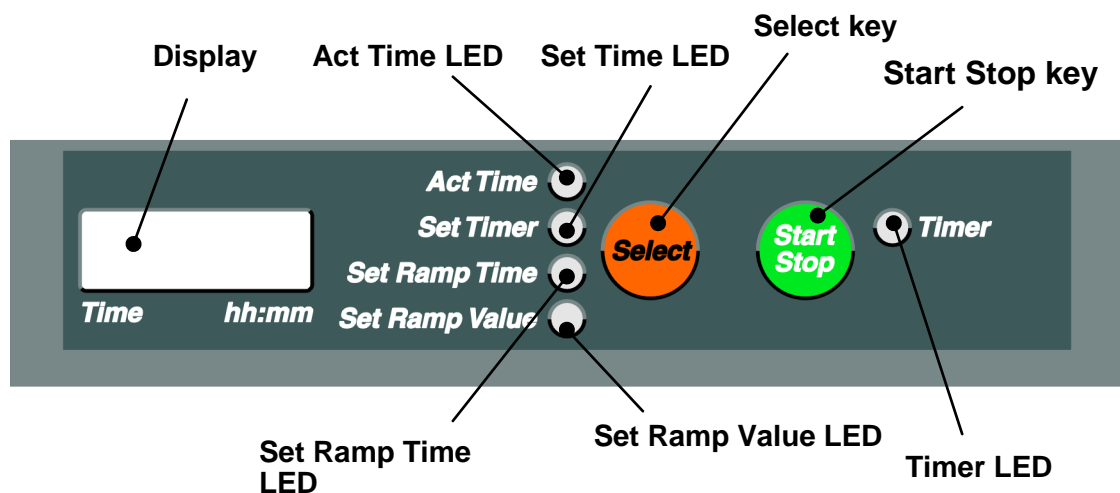

The timer function allows to end distillation after a preselected period of time has expired. After the time has elapsed, the following actions will be taken:

- Lift DOWN
- Rotation OFF
- Heater OFF
- Vacuum OFF
- Buzzer sounds for 5 seconds

increase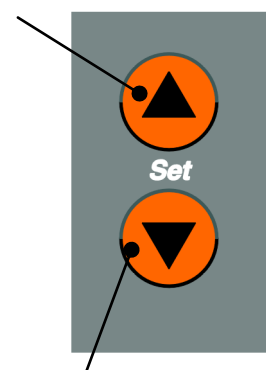

decrease

Turn timer ON and start

Select Timer with Select key.

Set hours and minutes with Set keys.

Start timer with Start Stop key, Timer LED flashes

Time remaining appears in display

Turn timer ON and start, after having started with Auto Start Stop before

Pressing **Start Stop** Timer, **Set Timer LED** illuminates.

For more settings refer to above paragraph.

19. Safety cut-out in an overpressure situation

The LABOROTA 20 control has been designed for distillation under vacuum. All of its gaskets and seals have been optimized for this vacuum atmosphere.

In particular situations, i.e. distillation in absence of a vacuum pump (e.g. low boiling substances), overpressure may build-up in the glassware kit. Please understand that our sealing system has not been designed for such overpressure situations.

In adverse overpressure situations, solvent may get into the drive module and hence damage bearings. As a safety precaution, when distilling in absence of vacuum, make sure to keep the aerating valve and valve (3) of the inlet pipe (2) open (also refer to "SET-UP", Chapter 8 page 74). Moreover LABOROTA 20 control features an electronic overpressure monitoring circuit; "Emergency function" will disconnect the item to avoid overpressure damage.

Description:

A vacuum sensor, installed to display and regulate distillation pressure, monitors pressure in the glassware kit either. Pressure rise above e.g. 1100 mbar will disconnect the LABOROTA 20 control by activating the „Emergency“ function:

Hot bath lift rises automatically,
 heater and rotation are disconnected,
 system is aerated,
 a buzzer sounds continuously,
Emergency LED lights, **Akt Vacuum LED** flashes.

For remedy choose distillation under vacuum or keep aerating valve and valve (3) on inlet pipe (2) open.

To reset Emergency-function, turn power switch OFF and ON.

Limit value:


Depending on local atmospheric pressure, limit value may be adjusted. In the factory, a standard value of 1100 mbar was set, you may do your own settings between 900 and 1200 mbar.

Rule of thumb:

The higher your local elevation, the lower your limit value. Make reference to the pressure in the aerated system, as displayed, plus about 50 mbar.

Edit limit value:

Turn LABOROTA 20 control OFF with "**power**" switch. Depress **Select** and **Start Stop** keys in the „Time“ block the same time, hold depressed and turn **power** switch to "ON". Wait for all of the **LEDs** are lighted and release **Select** and **Start Stop** keys.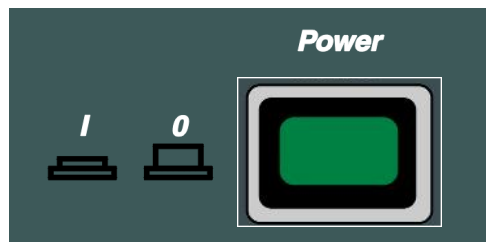

Select key

Start Stop key


„OVER“ appears in the **Display Speed**


E

Display I shows the actual limit value, e.g. 1,100 mbar. Use **arrow keys** to edit this value.

increase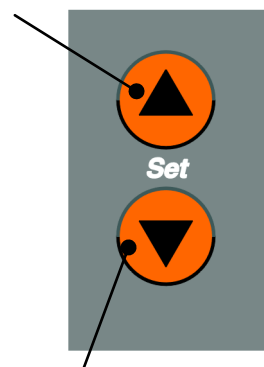

decrease

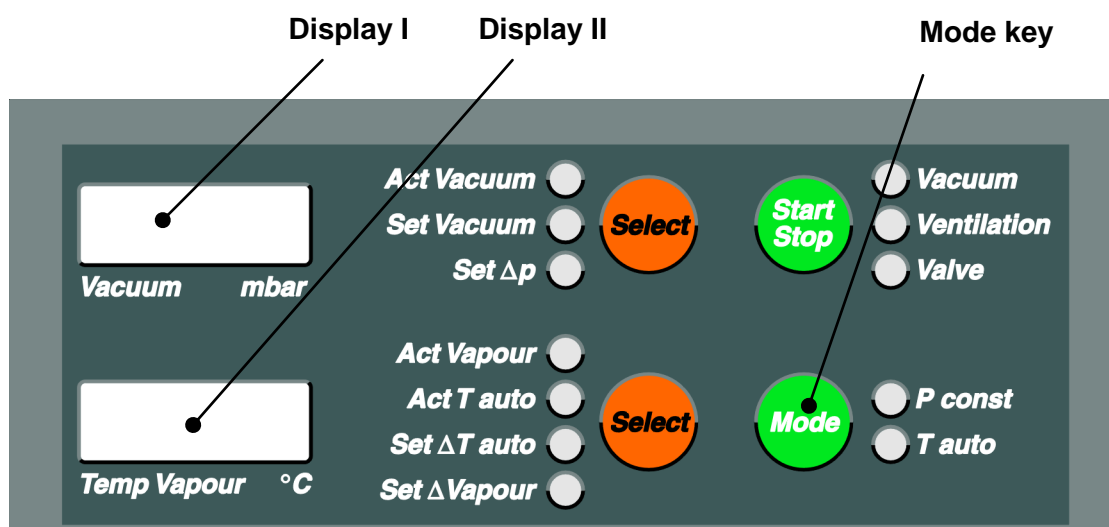Display II shows the emergency cut-out function active or not by „ON“ or „OFF“. Use **Mode** key to select „ON“ or „OFF“.

Do not override this safety feature except in very rare cases ! Your LABOROTA 20 might be damaged !

Confirm this action by turning the **power** switch OFF.

20. Safety cut-off in case of a defect in the vacuum system


The LABOROTA 20 control has a special monitoring function for the vacuum. Disconnected vacuum hoses, broken glass parts, defective vacuum pump, and lack of vacuum from the central vacuum supply are recognized during distillation. These are recognized when a previously set pressure limit value is exceeded. When this limit value is exceeded, distillation is stopped by the „Emergency“ function.

Description:

The vacuum sensor, which is installed to display and regulate the distillation pressure, is used for monitoring the pressure in the glass set. If the pressure rises above a limit value of 900 mbar, for example, with the vacuum regulation activated, distillation is stopped by the „Emergency“ function:
 The heating bath lift automatically moves downward,
 The heater and rotation stop,
 The system is aerated,
 Continuous buzzer,

Emergency LED illuminates, Akt Vacuum LED blinks, and “SAFE” is shown in the Vacuum Display alternating with the actual pressure.

Reset the Emergency-Function by switching the master switch on and off again.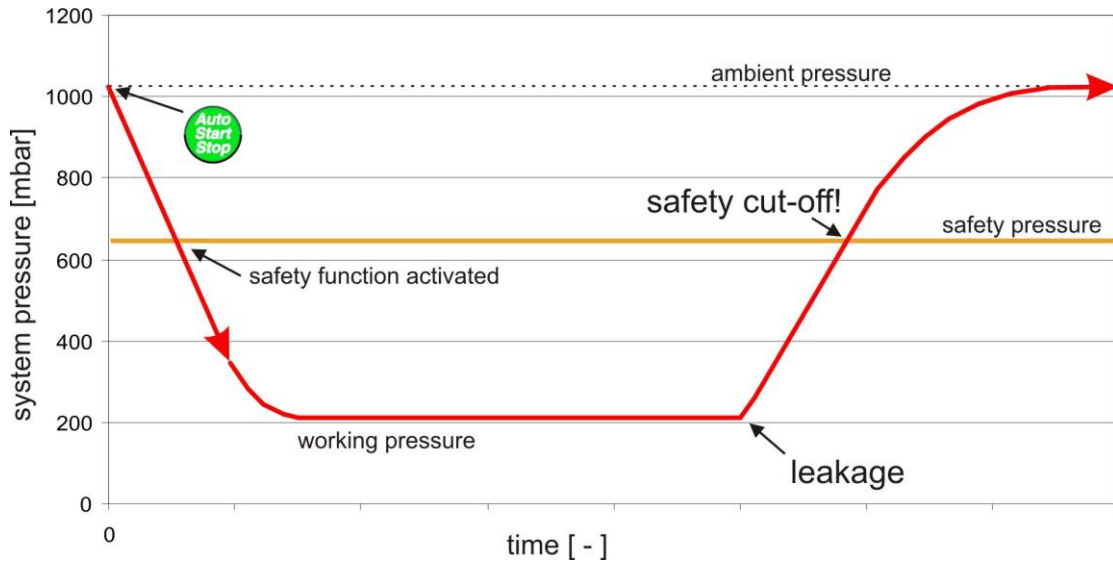

Limit value:

Depending upon the desired vacuum, at which distillation should occur, it can be advantageous to adjust the limit value. From the factory, the limit value is set at 900 mbar. The limit value can be changed from 0 to 1200 mbar.


Pay attention that the limit value is never set lower than the value, at which distillation should occur, as an „Emergency“ will be triggered otherwise.

Change limit value:


Switch off the LABOROTA 20 control with the master switch „Power“.

Simultaneously press the Select key and the Start Stop key in the „Speed“ field and switch on the master switch „Power“. Wait until all LED’s light up and then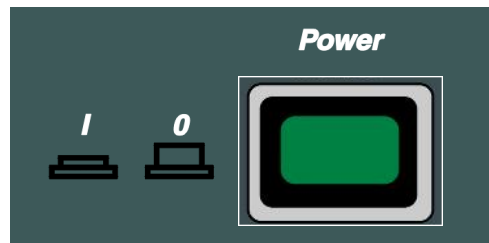

E

release the Select key and the Start Stop key.

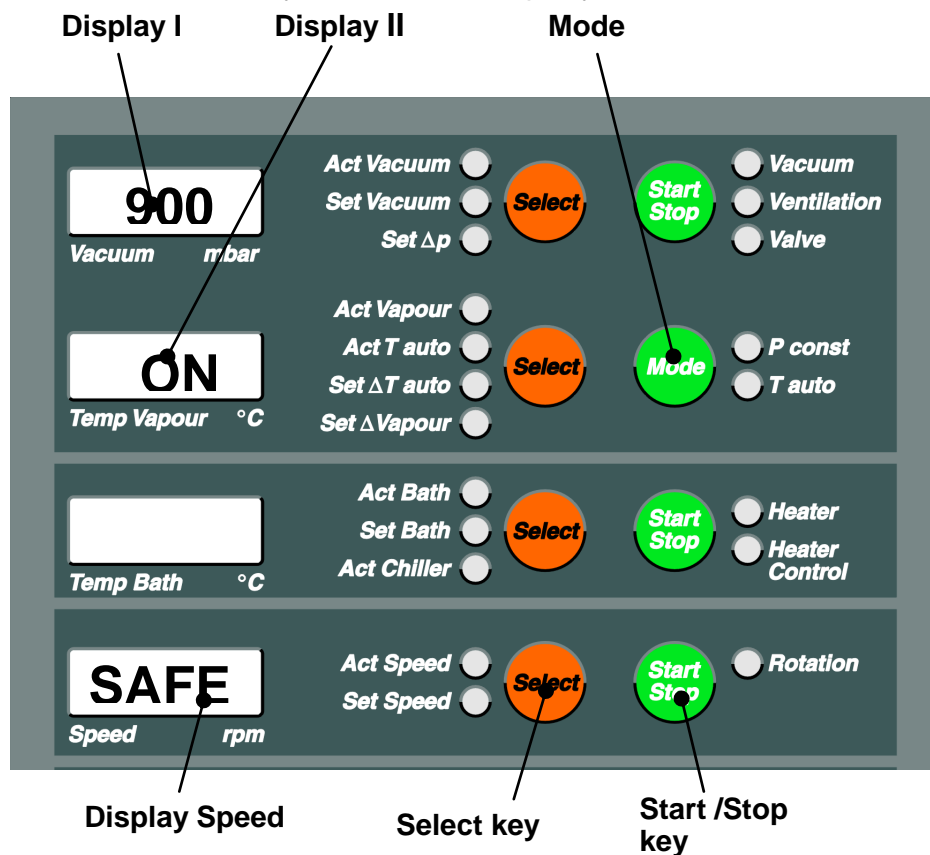„SAFE“ appears in the **Display Speed**

Whether or not the safety cut-off is active (ON) or switched off (OFF) is shown in **Display II** (Temp Vapour) with „ON“ or „OFF“. You can change from „ON“ to „OFF“ with the **Mode** key.

Switching off the safety cut-off is the sole responsibility of the user.

This function does not work during distillation without vacuum! Here, the safety cut-off offers protection in case of excess pressure.


Even when using the auto function, the function only activates after the limit value is exceeded. Ensure by experimentation that the set limit values are never reached!

The current limit value is shown in **Display I** (vacuum) e.g. 900 mbar. The limit value can be changed with the **arrow** keys.

The process must be concluded by switching off the master switch „**Power**“.

Increase value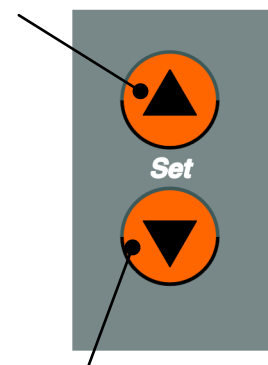

Decrease value

21. Safety functions

All of the following cause LED ON situations.


- The **Safety Cover LED** tells you that the enclosure is open. Rotation won't start, or rotation is going to stop. Close hood and hit Start Rotation key.
- **Flask Support LED**: Flask support is not in its extreme position (either side). Rotation won't start, or rotation is going to stop, lift won't work. Move flask support to extreme position (both sides) and secure with knurled knobs (5), (page 103).
- **LED Rotation Stop LED**: Drive lock engaged, rotation can't start. Release lock by turning knob CCW.
- **Overheat Protection LED**: Bath temperature beyond acceptable limits. Heater won't start, heater is disconnected. Check heater for possible problems.
Overheat Protection LED flashing: After response, over temperature protection enters Emergency mode (refer to next paragraph).
 Heater won't start, heater is disconnected.
 Possible reasons: heating bath level insufficient, heating bath dry; refill.
 Heater is broken.
 After repair action allow bath to cool off and reset over temperature protection by depressing little button with screwdriver.
- **Emergency LED**: Emergency CUT-OUT actuated or over temperature protection response (refer to para. above)
 Heating bath lift descends automatically;
 Heater and rotation OFF;
 vacuum is turned OFF and system aerated.
 Continuous buzzer
 Reset with master switch OFF and ON
- The **Emergency LED** being lighted and the **Akt Vacuum LED** flashing the same time means the safety circuit detected an overpressure situation (refer to para. 19 page 90).
 Heating bath lift descends automatically;
 Heater and rotation OFF;
 vacuum is turned OFF and system aerated.
 Continuous buzzer
 Reset with master switch OFF and ON

TIMER AND RAMP-FUNCTION

Timer and ramp functions allow to program time ramps for vacuum pressure (p const mode), ΔT auto (T auto mode), bath temperature and speed. 10 steps are available for every parameter. A total of 40 individual times can be programmed.

22. Set ramp values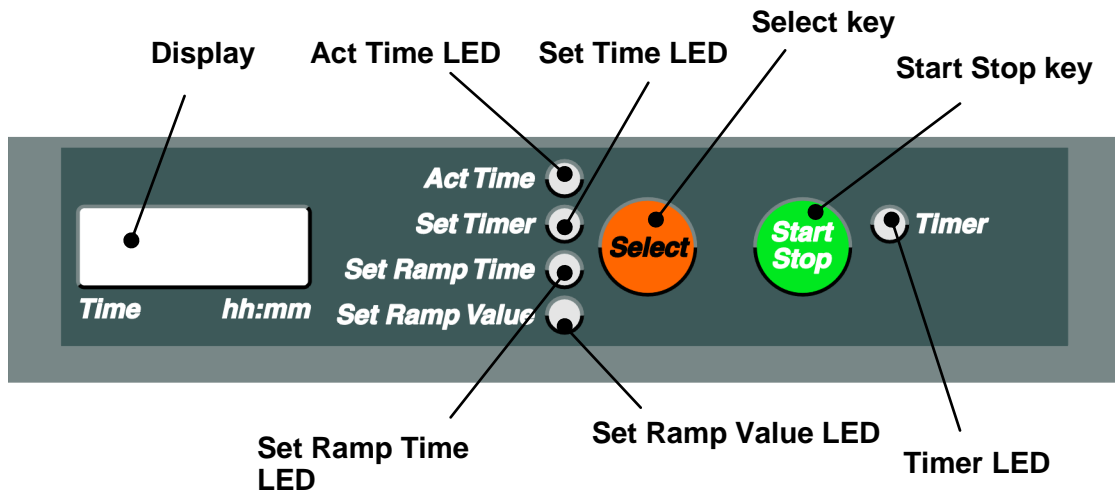

Before programming a new ramp, make sure to have deleted the old one (refer to “delete ramp”, page 100).

Procedure 1st step:

You need to program 00:00 for the first time.

Press the **Select key** next to the **Display Time** until **Set Ramp Time LED** illuminates.

Set Ramp Time LED ON. **Display Time** only is active, all other displays are blank.

Display shows time for the present ramp in hh:mm.

The time 00:00 should always be programmed as the first time (is preset).

Press the **Set keys** to alter time setting (**forward & back**).

Not pressing a key for 5 seconds, display will return to Act. Last value entered will be saved.

To program parameter values as a function of times selected, go to **Set Ramp Value LED** with the **Select key** next to the **Display Time**. **Set Ramp Value LED** illuminates, **Set Ramp Time LED** flashes.

Go to parameter desired with **Select key** next to the **Display Time**, **Set LED** illuminates: the following parameters can be programmed:

vacuum (p const mode) **Set Vacuum LED** illuminates

ΔT auto (T auto mode) **Set ΔT auto LED** illuminates

bath temperature **Set Bath LED** illuminates

speed **Set Speed LED** illuminates

If one or more of the parameters listed above is not to be assigned a ramp function, leave this parameter as „OFF“.


As shipped, no ramp functions are set. (“OFF” displayed)

Values are programmed with the **Set keys**, as described.


Write value in memory by pressing the **Select key** once more.

Use the same procedure to program the next parameter.

increase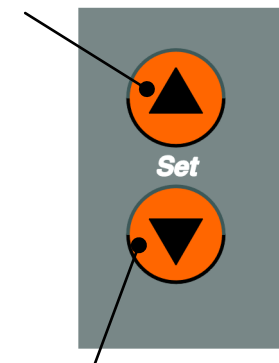

decrease

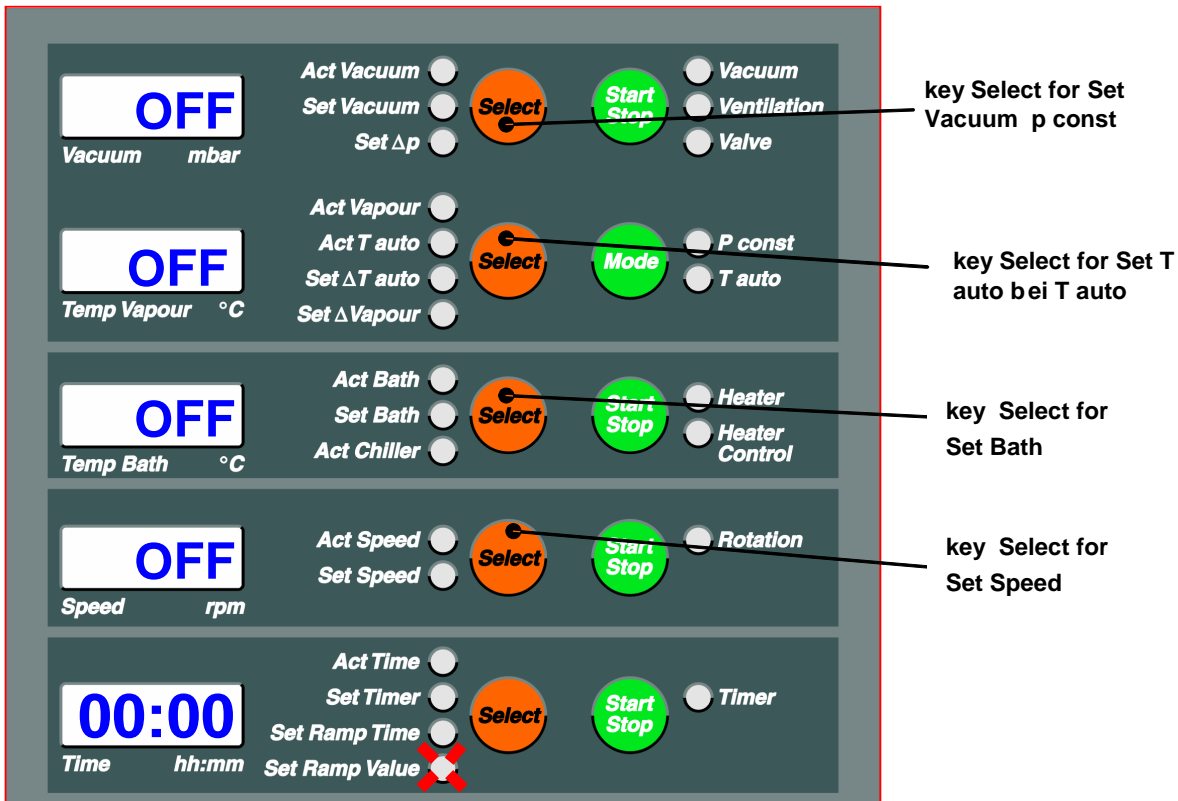2nd step:

Now, select the LED Set Ramp Time by pressing the **Select key** next to Display Time repeatedly. The time for the 2nd step of programming can be changed with the Set increase or Set decrease keys.

After 5 seconds with no entry, the display returns to Act Condition. The last value is accepted.

Additional steps:

Additional steps can be programmed in the same manner as step 2.

Select **Set Ramp Time LED** to program the next time.

For complex ramps, we recommend you to use the form enclosed (for convenience). The same page shows an example.

Note: Values only can be programmed that had been activated at 00:00 before.

Parameters, which were programmed as a ramp, cannot be changed during the process of the ramp with the „Set Function“. Parameters that are not programmed as a ramp can be changed at any time with the „Set Function“.


Time	Unit	1	2	3	4	5	6	7	8	9	10
Set Ramp Time	hh:mm										
Set Vac	mbar										
Set T auto	°C										
Set Bath	°C										
Set Speed	rpm										

Time	Unit	11	12	13	14	15	16	17	18	19	20
Set Ramp Time	hh:mm										
Set Vac	mbar										
Set T auto	°C										
Set Bath	°C										
Set Speed	rpm										

Time	Unit	21	22	23	24	25	26	27	28	29	30
Set Ramp Time	hh:mm										
Set Vac	mbar										
Set T auto	°C										
Set Bath	°C										
Set Speed	rpm										

Time	Unit	31	32	33	34	35	36	37	38	39	40
Set Ramp Time	hh:mm										
Set Vac	mbar										
Set T auto	°C										
Set Bath	°C										
Set Speed	rpm										

Time	Unit	1	2	3	4	5	6	7	8	9	10
Set Ramp Time	hh:mm	0:00	0:15	0:20	0:30	0:40	0:50	1:00	2:00		
Set Vac	mbar	1000	1000	700	400	200	150	100			
Set T auto	°C	OFF									
Set Bath	°C	20	50								
Set Speed	rpm	20	30								

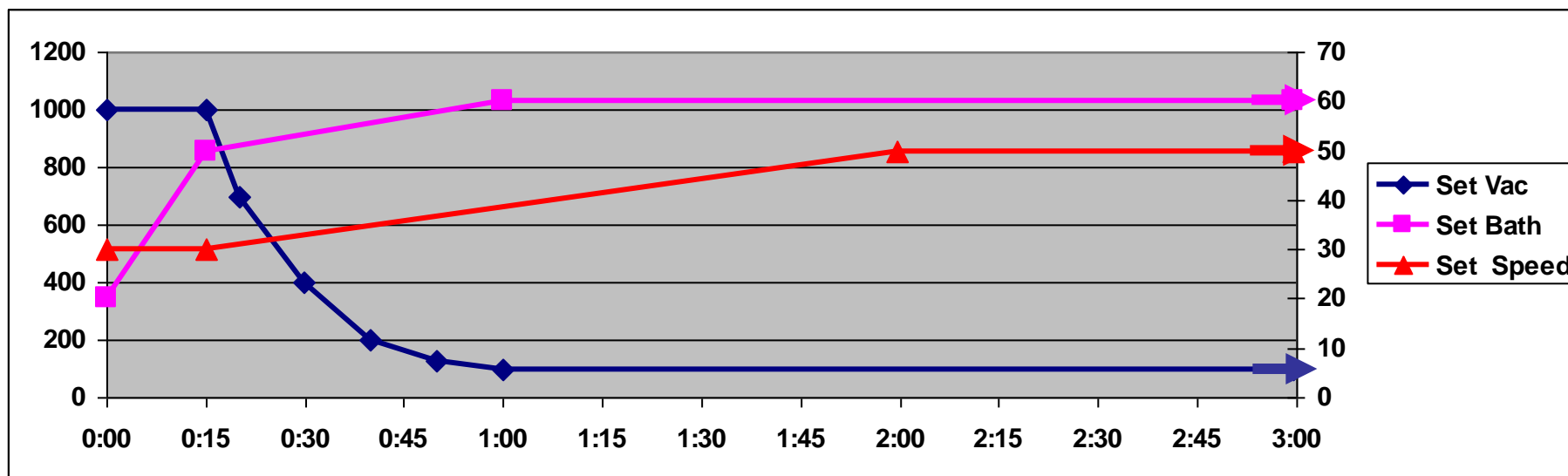23. Display ramp values

Ramp values once programmed can be displayed and edited.

Hit **Ramp key**, **Ramp on LED** illuminates. Press the **Select key** (Time) until **Set Ramp Value LED** illuminates.

First programmed time 00:00 and corresponding value are displayed.

Void parameters are marked OFF.

More times and programmed values are displayed with the Set keys.

24. Edit ramp values

To edit a value, select with **Select key**.

Now edit value with the **Set keys**.

Accept new value by pressing the **Select key** once more.

Program more times: refer to para. 1, set ramp values.

25. Start ramp function

Program ramp values as described in 1).

Press **Ramp key**, **Ramp on LED** illuminates.

Hit **Auto Start Stop key**, **Auto Start Stop LED** illuminates.

The following functions are started:

lift UP

rotation ON

vacuum ON

heater ON

All values programmed for a ramp are controlled automatically.

Display shows time elapsed from START.

Values between times programmed are interpolated.

26. Discontinue ramp function

Ramp function may be discontinued any time.

To discontinue hit the **Ramp key**, **Ramp LED** dims.

The very last ramp function values are kept.

A ramp once programmed will be kept in memory even during power-off times and can be used and edited all the time.

27. Delete ramp

Ramp LED and Set Ramp Value LED must illuminate.

Press both **SET keys** at the same time and hold. After 5 seconds, "del" appears for 3 seconds, a buzzer sounds.

The ramp has been deleted.

With the ramp function activated, it is not possible to also activate the timer function! If the timer function is activated during the ramp function, the ramp function will be switched off and the current values of Set Vac, Set Bath, Set Speed, and Set ΔT auto are accepted as Set-values. These can then naturally be changed.

INTERFACE

28. Connect interface

Make sure the Unit is OFF before you attempt to connect the interface cable.
Use a cable with SUB-D 9 connectors and connect with the terminal connector located underneath the **LABOROTA 20 control** control panel.

29. Interface parameters

9600 Baud, no parity, 8 bits, 1 stop bit, no protocol

30. RS 232 interface

This interface allows you to specify all Set-values, move the lift up and down, open and close aerating valve and vacuum valve.

All Act-values and status of safety functions may be recalled.

31. RS 232 interface commands

query

Command	action	remarks
Act commands		
VAC?	Act Vacuum	
VAP?	Act Vapor	
TAU?	Act T auto	
BATH?	Act Bath	
CHILL?	Act Chiller	
SPEED?	Act Speed	
TIME?	Act Time	elapsed time, continuous operation
COUNT?	Act Time	time, timer mode
SET1?		query Set-value
SET2?		query Act-value
Safety functions		
SC?	Safety Cover	0 = action; 1 = no action
FS?	Flask Support	0 = action; 1 = no action
RS?	Rotation Stop	0 = action; 1 = no action
OP?	Over Temperature Prot.	0 = action; 1 = no action

Triggering

Command	action	remarks
Set commands		
VAC=XXXX!	Set Vacuum	4-digit in mbar; range set 0000 to 1200
DP=XXX!	Set Δp	3-digit in mbar; range set 000 to 200
DT=XXX!	Set ΔT auto	3-digit, one fractional part, in °C range set 001 to 300 equals 0.1 to 30.0
DV=XXX!	Set Δ Vapor	3-digit, one fractional part, in °C range set 001 to 200 equals 0.1 to 20.0
BATH=XXXX!	Set Bath	4-digit, in 1°C increments range set 0000 to 1800 equals 0.0 to 180.0
SPEED=XXX!	Set Speed	3-digit, in rpm; range set 000 to 180
Actions		
SVX!	Start Stop Vacuum	SV0! = OFF ; SV1! = ON
VVX!	Valve	vacuum valve; VV0! = CLOSE ; VV1! = OPEN
BVX!	Ventilation	aerating valve; BV0! = CLOSE ; BV1! = OPEN
MP!	Mode p const	go to p constant mode
MT!	Mode t auto	go to T auto mode
SBX!	Start Stop Bath	SB0! = heater OFF ; SB1! = heater ON
SSX!	Start Stop Speed	SS0! = rotation OFF ; SS1! = rotation ON
LU!		lift UP
LD!		lift DWN

FUNCTIONS AND HANDLING


32. Protective enclosure

The protective enclosure (1) is an extra safety feature for the operator. Its large pane of a special safety glass guarantees a clear view of the rotating flask. A handle (2) on the front is used for closing and opening. In both extreme positions, it is held by a pneumatic strut.

Caution: Flask will not rotate unless the hood is closed. Opening the enclosure during operation will stop rotation.

Safety Cover LED (6) illuminates.

After closing the enclosure, rotation is restarted by the Start Stop key.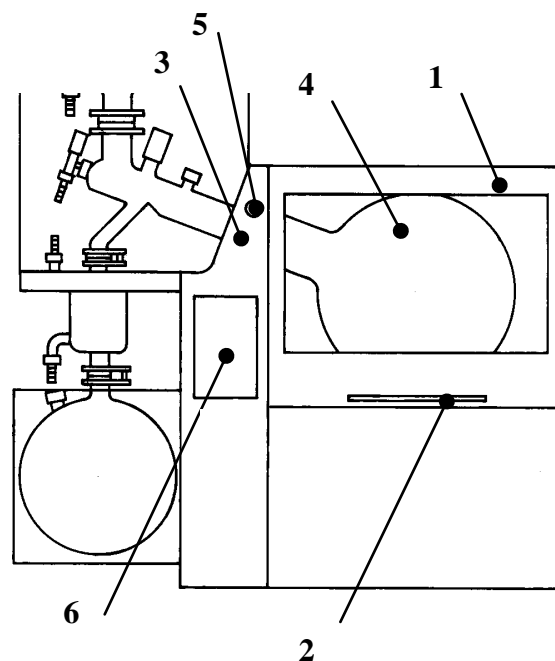

33. Drive lock

picture on page 89

The drive lock facilitates opening and closing the large coupling ring (3), holding the flask (4).


- To engage lock: turn knob (5) CW until “**lock**” appears.
Manually rotate drive head, until lock snaps into place. **Rotation Stop LED** (6) on front panel illuminates.
- To release lock: turn knob (5) CCW until “**unlock**” appears.
Rotation Stop LED (6) on front panel dims.

34. Tightening of Evaporator Flask


Evaporator Flask must be sufficiently tightened with large coupling ring (X).

For this purpose the large coupling ring (X) is equipped with a torque indicator. As illustrated below the coupling ring has to be tightened until the indicator point (Y) is aligned with point „1“ (Z). This is accompanied by a „click“. For distillation without vacuum the ring has to be tightened until point (Y) is aligned with point „2“ (A).

If rule is not followed leakage can occur during operation.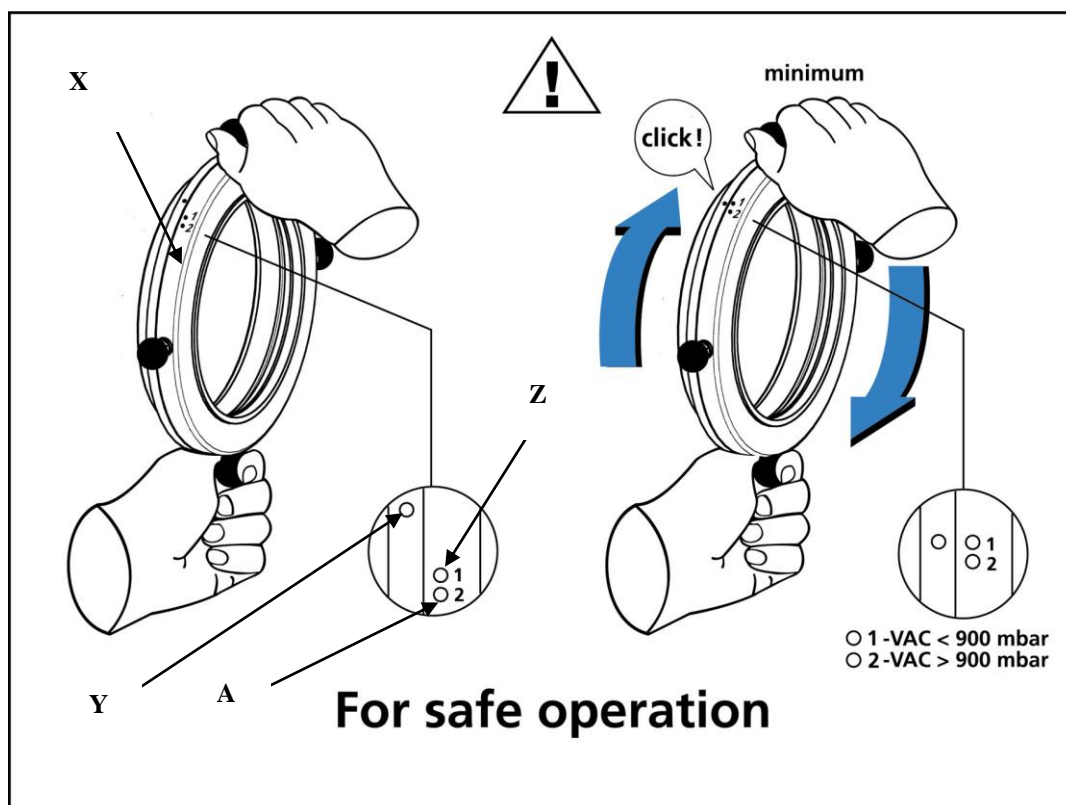

35. Flask supports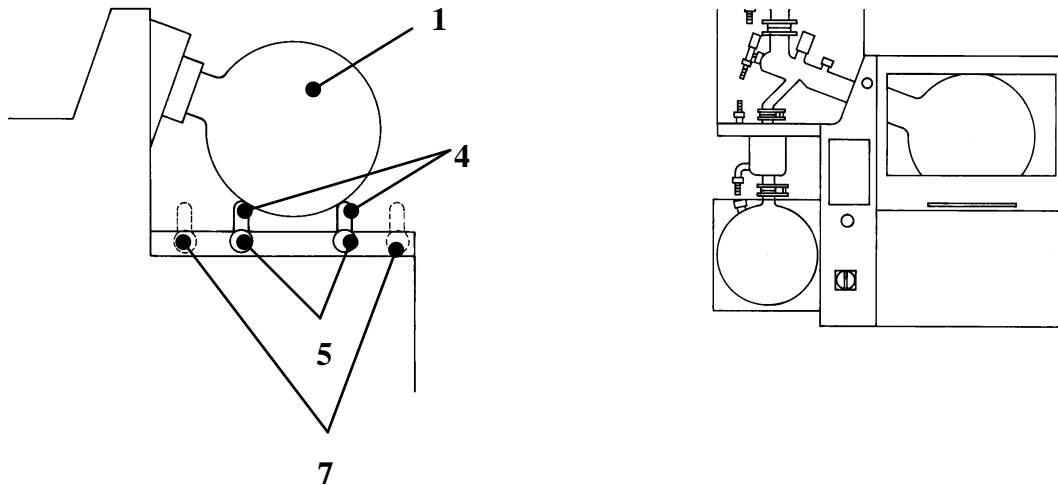


The flask supports make flask (1) removal (even filled) an easy job for one person.

3. Uninstall flask: Open enclosure (2). Engage drive lock (3). Slide support pipes (4) to flask (1) from both sides. Check for firm contact and lock support pipes in place by tightening thumbscrews (5) CW. Unscrew large coupling ring (6, not shown) by turning CCW. Tilt flask neck with large coupling ring (6) forward. Remove flask carefully (coupling ring to remain on flask).
4. Install flask: Open enclosure (2). Engage drive lock (3). Move support pipes (4) (both sides) to the respective flask size mark (center of thumbscrew = mark). Mark I = 10 l flask; mark II = 20 l flask. Lock support pipes in place by tightening thumbscrews (5) CW. Put flask (1) , holding large coupling ring (6), on support pipes carefully. Swing flask axis with coupling ring in-line with axis of rotation, until flange is centered in drive head. Secure PTFE-adapter (not shown) with PTFE-O-ring tight by CW turn. Double-check for firm seat. Return support pipes to initial position (7). Release drive lock (3). **Caution:** Basic position of support pipes is monitored by the electronic circuit; when leaving this position, all electric functions of the system will freeze. This situation is reported by the Flask Support LED on the control panel.

36. Automatic water refill (water bath model only)

The LABOROTA 20 control WB features a automatic water refill system with float switch.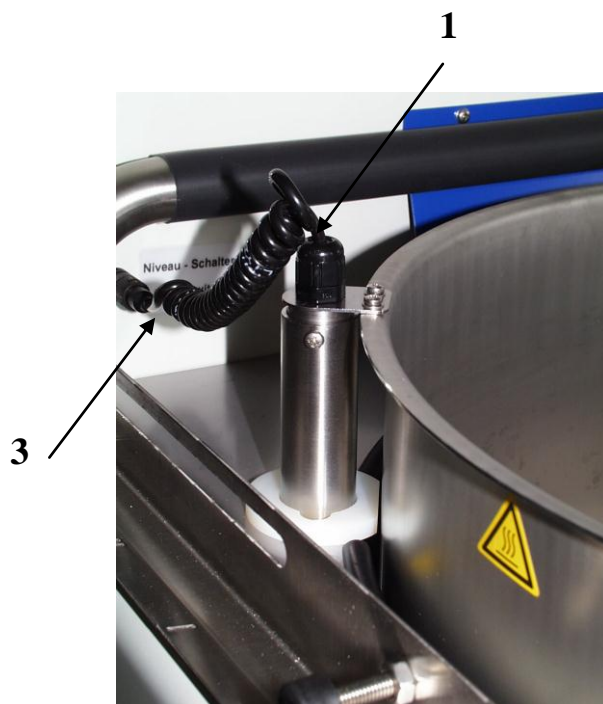

The float switch (1) monitors water level in the heating bath. When dropping below normal level (about. 30 mm below tank rim), a solenoid valve (12, next para.) opens the water supply until reaching the normal fill level. Cooling water is supplied to the heating bath via pipe. The float switch needs to be connected with plug connector (3).

Caution: For technical reasons automatic refill requires the heating bath lift to be at its upper most position.

Upward lift travel is limited by the float switch. Water control valve on the rear must be connected.



The float switch may fail due to calcareous deposits, causing the hot bath to be overfilled or overflow. We recommend to decalcify this float switch every 6 months.

To this end uninstall float switch (disconnect plug connector (3) and loosen float switch (1) attaching screws with a screwdriver.

Use commercial grade decalcifying agent (go by instructions on the pack), and immerse float switch until fluid level reaches coupling nut. For treatment time refer to instructions on the pack.

37. Vacuum and water distribution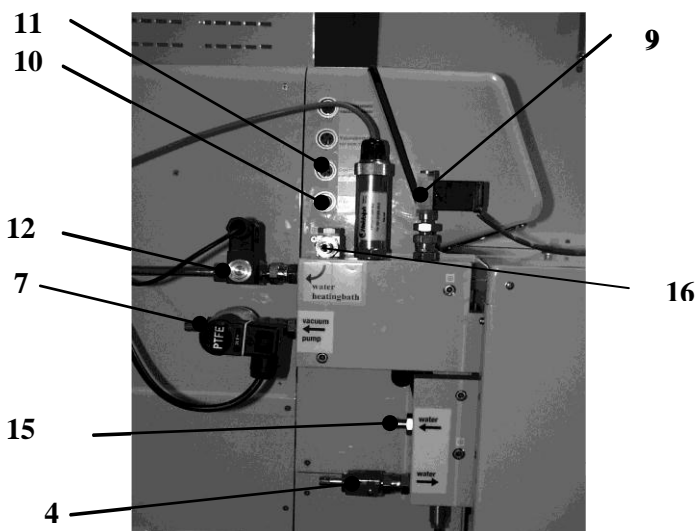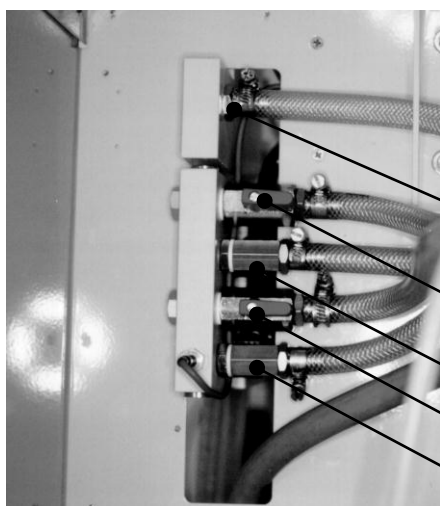

The rear panel of the unit contains a vacuum/water distributor. A central cooling water supply optionally works with:

1 chiller ; 1 chiller plus 1 after-chiller, or 2 chillers. Refer to picture when connecting the water hoses (ID=8 mm). Hoses are routed in the chiller panel and attached to the chiller with hose nipples. Inlet flow control uses a ball valve (4). When using 2 chillers, open ball valve (5) (lever in-line); when using 1 only chiller, close ball valve (6) (lever out-of-line). The same applies to the coolant return network. When using 2 chillers, open ball valve (6) (lever in-line), using 1 only chiller, close ball valve (6) (lever out-of-line).

Connect #1 chiller with the fittings (13). Coolant return is connected with fitting (15).

Vacuum pump is connected with vacuum fitting (7) (shown with optional vacuum valve), use ID = 8 mm vacuum hose). Connect vacuum distributor with fitting (8) on glass kit. Vacuum hose is routed in the chiller panel to glass kit fitting and attached with screwed nipples. The built-in aerating valve (9) is controlled from the control panel.

The aerating valve (9) is connected with #3 terminal connector (6-pin) (11).

Water refill solenoid valve (12) is connected with the lower terminal connector (4-pin) (10).

The cooling coils of the cooler may only be exposed to a maximum cooling water pressure of 2 bar. So that higher pressure can never develop, no valve and no restrictor may be installed in the cooling water return.

Always secure the cooling water hoses with hose clamps!

38. Pt100 temperature sensor, chiller

For T auto mode, a chiller Pt100 temperature sensor needs to be connected with terminal connector (1).

HEATING BATH FUNCTION AND HANDLING

39. Fill heating bath (water bath only)

Using de-ionized or distilled water needs to add 0.2 % of Borax ($\text{Na}_2\text{B}_4\text{O}_7 \cdot 10\text{H}_2\text{O}$) (corrosion inhibitor)

Please note: The use of tapwater as heating media can lead to calcification which can provoke crevice corrosion. Therefore the user is strongly recommended to clean the heating bath with an adequate polish on a regular basis.

Heating bath water supply uses the vacuum/water distributor (refer to page 105).


Connect inlet end to fitting. Heating bath fill uses line (12) and is controlled by a solenoid valve (12). Also refer to page 105.

40. Drain heating bath (water bath only)

The rear panel of the LABOROTA 20 control with water bath features a drain valve (1). To close valve turn black lever perpendicular to line.

This valve drains bath fill into a tank (not supplied).


Caution: Allow water to cool to ambient to avoid burns.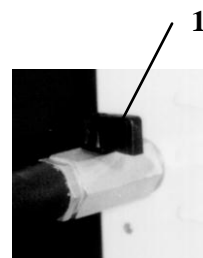

41. Condensate drain

The LABOROTA 20 control (water bath) features a condensate drain. Water vapor from the hot bath is condensed by the chiller coil (1) of the condenser (2), and is collected in a sump (3), which in turn is drained through a nipple (4).

This nipple (4) accepts an elastic hose (ID = 8 mm). The other end is plugged in a collection tank.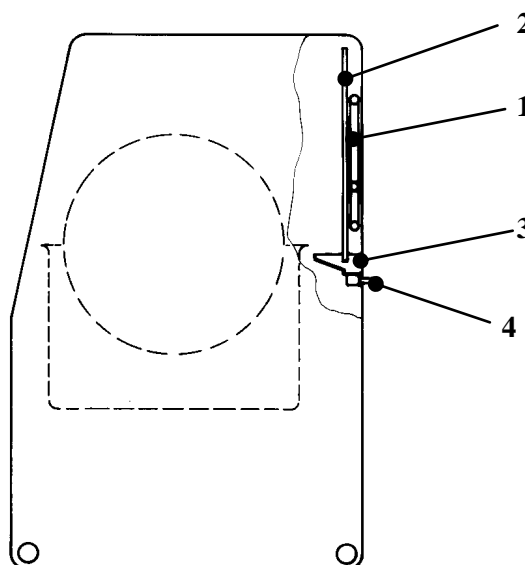

42. Heating bath over temperature protection

Once the over temperature protector responds (Over temperature Protection LED), allow the bath to cool off and fix the problem.

Most possible reason: no water supply.

Descend heating bath lift to LOW.


Turn master switch OFF.

Reset over temperature protector by introducing an insulated screwdriver through the little hole at the lower right end of the unit. (remove black plastic plug to gain access through panel)

43. Vacuum pressure sensor/vacuum valve


In order to be able to regulate the vacuum, the vacuum pressure sensor (1) (No.: 591 21000 00) and the vacuum valve (2) (No.: 591 20000 00) are already installed.

The vacuum pressure sensor is connected to the top terminal connector (3) and the vacuum valve to the second terminal connector (4).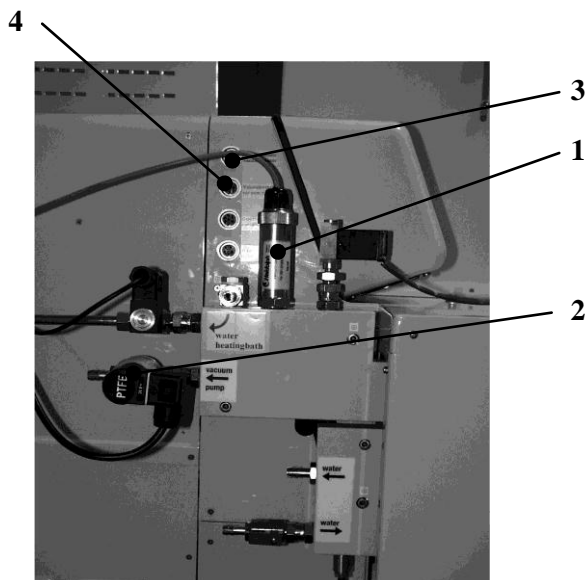

44. Temperature sensor for vapor


To display vapor temperature on the control panel, you'll need the optional temperature sensor (1) (P/N 14-014-003-06). To install, remove cap from temperature fitting (2) of glass kit. Its plug fits the rear #3 terminal connector of the drive head. Secure in place.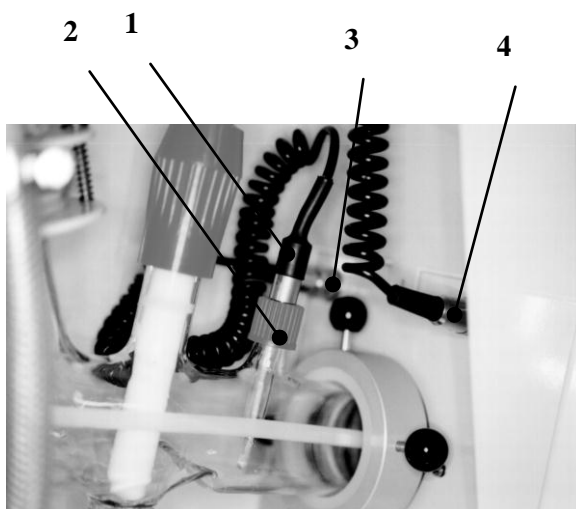

45. Temperature sensor for T auto vacuum mode

T auto vacuum control mode needs the optional temperature sensor (5) (P/N 14-014-003-06).

To install, remove caps from condenser (6). When running the T auto mode, connect temperature sensor with the forward terminal connector (4) in the rear of the drive head.

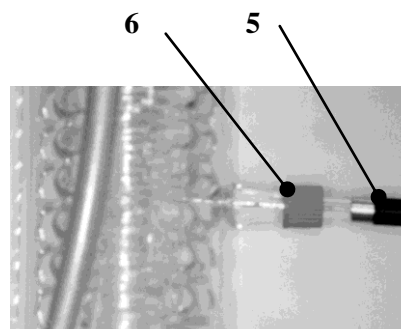CLEANING & SERVICING

Cleaning: wipe housing clean with a damp cloth (add some mild liquid soap).

Note

To avoid damage to the surface finish, avoid using chlorine bleach, chlorine-based detergents, abrasive substances, ammonia, rags or cleaning agents containing metal particles.

Rust stains appearing on the bottom of the hot bath after long-term use are caused by foreign matter deposits (iron particles or rust particles). Remove such rust stains and calcareous deposits with a conventional phosphoric acid bathroom cleaner (decalcifying agent).

The float switch may fail due to calcareous deposits, causing the hot bath to be overfilled or overflow. We recommend to decalcify this float switch every 6 months.

To this end uninstall float switch (disconnect plug connector (3) and loosen float switch (1) attaching screws with a screwdriver. Use commercial grade decalcifying agent (go by instructions on the pack), and immerse float switch until fluid level reaches coupling nut. For treatment time refer to instructions on the pack.

Never attempt to work with damaged or worn-out PTFE gasket (6). Not reaching vacuum pressure wanted may damage your rotary evaporator. Inspect PTFE gasket (6) in regular intervals and replace on condition.

The unit is maintenance-free. Repair work is limited to technicians approved or appointed by Heidolph Instruments. Please call your local Heidolph Instruments Dealer or a Heidolph Instruments Field Representative (also refer to page 111).

UNINSTALL, FORWARD & STORE

Uninstall

Turn the item OFF and disconnect mains plug.

Forward & Store

1. We recommend to store the item and its components in its original box, or a similar container that offers adequate protection against damage in transit. Tape the box securely.
2. Store the item in a dry place.

Caution

Do not jolt or shake the item during transport.

DISPOSAL

For disposal, please comply with your local or national regulations. Separate by metal, plastic, etc. Packing material to be treated as described above (material split).

TROUBLESHOOTING

- Master switch won't light when turning ON
check power cord
master switch broken (call service)
- Display: Act Vacuum "0000"
vacuum sensor broken or not connected
- Display: Act Vacuum won't change during evacuation
vacuum valve connected with vacuum pressure sensor terminal
- Display: Act Vapor "----"
temperature sensor not connected or broken
- Display: Act T auto "----"
T auto temperature sensor not connected or broken
- Display: Act Chiller "----"
Cooling water temperature sensor not connected or broken
- Vacuum valve inoperative
vacuum pressure sensor not connected with valve terminal, or valve broken
- Display: Act Vacuum (aerated) differs from atmospheric
altitude correction factor not integrated.
Calibrate vacuum pressure sensor.
- Poor vacuum pressure
inspect gaskets and fittings
check ground fittings, coat with vacuum grease, as required
PTFE seal installed the wrong way. Sealing lip needs to point towards drive head
check line to vacuum pump
- Heating bath lift won't travel up
float switch not connected or broken
also refer to safety functions section
lift mechanism broken
upper limit switch broken
- Heating bath lift won't travel down
lift mechanism broken
lower limit switch broken
- No rotation
also refer to safety functions section
motor broken
- Mode won't change from P const to T auto.
T auto temperature sensor not connected or broken
- Aerating valve won't work. Aerating valve broken or not connected.
- **Distillation automatically discontinued (by accident)**
EndP value incorrect; **Set Time** incorrect, cut-out sensor responded.

Also refer to safety functions (page 95)

Whenever facing a problem that cannot be fixed by the above information, call your authorized Heidolph Instruments dealer right away.

WARRANTY, LIABILITY & COPYRIGHT

Warranty

Heidolph Instruments provides a three-year warranty on the products described here (with the exception of glass and consumable parts) if registered with enclosed warranty card or via internet (www.heidolph.com). Warranty starts with the date of registration. Without registration warranty starts according to serial number. This warranty covers defects in materials and workmanship. Transit damage is excluded from this warranty.

To obtain such warranty service, contact Heidolph Instruments (phone: +49 – 9122 - 9920-68) or your local Heidolph Instruments Dealer. If defects in material or workmanship are found, your item will be repaired or replaced at no charge.

Misuse, abuse, neglect or improper installation are not covered by this warranty promise.

Alterations to the present warranty promise need Heidolph Instruments' consent in writing.

Exclusion Clause

Heidolph Instruments cannot be held liable for damage from improper use or misuse. Remedy for consequential damage is excluded.

Copyright

Copyright in pictures and wording of the present Instruction Manual is held by Heidolph Instruments.

QUESTIONS / REPAIR WORK

If any **aspect** of installation, operation or maintenance remains unanswered in the present Manual, please get in touch with the following address:

For repairs please call Heidolph Instruments (phone: ++49-9122-9920-68) or your local, authorized Heidolph Instruments Dealer.

Note

You will receive approval for sending your defective item to the following address:

Heidolph Instruments GmbH & Co. KG
Lab Equipment Sales
Walpersdorfer Str. 12
D-91126 Schwabach / Germany
phone: ++49-9122-9920-68 Fax: ++49-9122-9920-65
E-Mail: Heidolph.Instruments@Heidolph.de

**Safety Information**

When shipping items for repair that may have been contaminated by hazardous substances, please:

- advise exact *substance*
- take adequate *protective* action towards our parts receiving and service personnel
- *mark* the pack IAW Hazardous Materials Act

**CE-DECLARATION OF CONFORMITY**

We herewith declare that the present product complies with the following standards and harmonized documents:

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© Heidolph Instruments GmbH & Co. KG